





Cellvento® and EX-CELL® Feed Mixing Protocol

1.

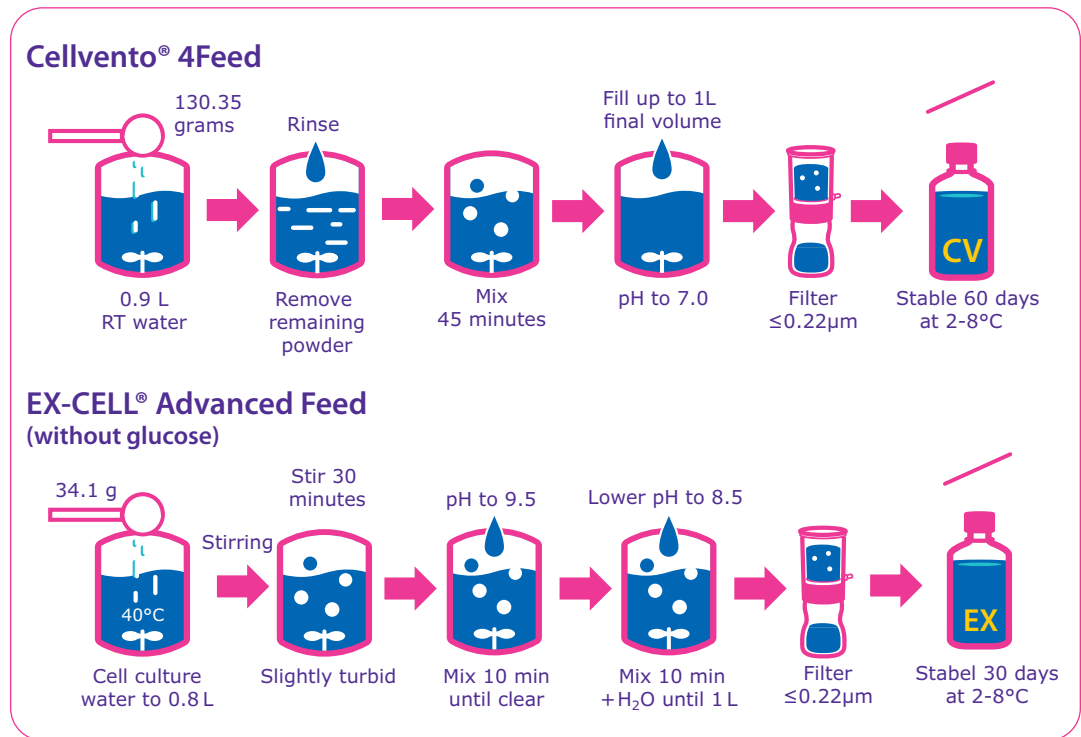

2.

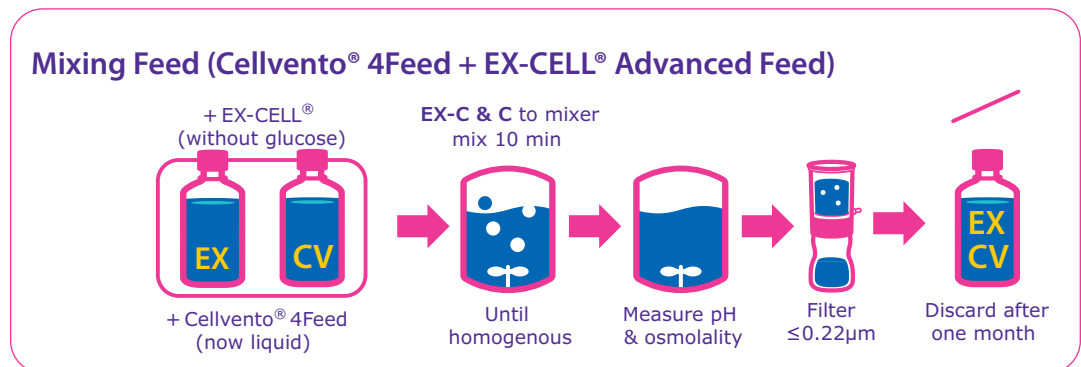

Additional information on back side.

Try different mix levels to find the optimal level of feed for your process

Suggested Mixing Levels

Mix Ratio	Cellvento® 4Feed	EX-CELL® Advanced Feed
Volume for one liter of mixed media		
100	1 L	0 L
75:25	750 mL	250 mL
50:50	500 mL	500 mL
25:75	250 mL	750 mL

Evaluate Cellvento® 4Feed as a mix and a standalone feed. Usually EX-CELL® Advanced Feed performs best with the Cellvento® 4Feed.

In some applications, Cellvento® 4Feed may perform better without EX-CELL® Advanced Feed.

To place an order or receive technical assistance

In Europe, please call Customer Service:

France: 0825 045 645

Germany: 069 86798021

Italy: 848 845 645

Spain: 901 516 645 Option 1

Switzerland: 0848 645 645

United Kingdom: 0870 900 4645

For other countries across Europe, please call: +44 (0) 115 943 0840

Or visit: [MerckMillipore.com/offices](https://www.MerckMillipore.com/offices)

For Technical Service visit:
[MerckMillipore.com/techservice](https://www.MerckMillipore.com/techservice)

[MerckMillipore.com](https://www.MerckMillipore.com)

Merck KGaA
Frankfurter Strasse 250
64293 Darmstadt, Germany

