


# Buchwald G6 Precatalysts: Oxidative Addition Complexes for Efficient L-Pd(0) Generation





Buchwald G6 precatalysts and other OACs have been applied as effective catalysts for the formation of C-C, C-N, C-O, C-F, and C-S bonds.<sup>1-5</sup> For the following reactions and applications, see **Table 2** for referenced Buchwald G6 and other OAC precatalysts.

**Table 1.** Reactions and Applications










<b>Fluorination of Aryl Bromides<sup>1</sup></b>			
	<p>70%</p>	<p>90%</p>	<p>92%</p>
<b>Fluorination of Aryl Triflates<sup>1</sup></b>			
	<p>82%</p>	<p>87%</p>	<p>94%</p>

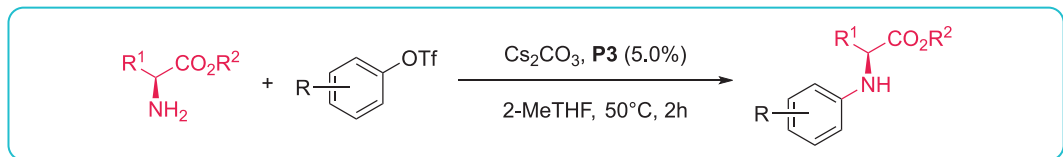






**Amino Acid Ester Arylation<sup>1</sup>**

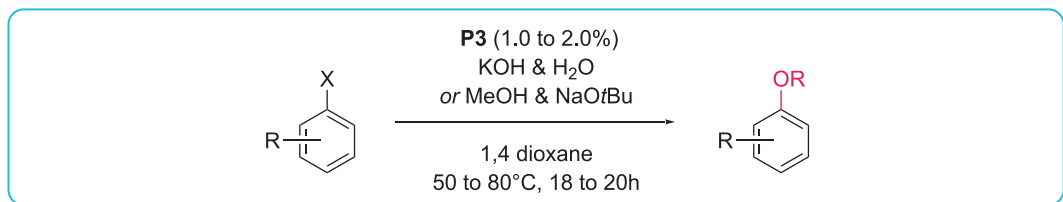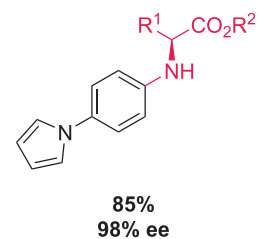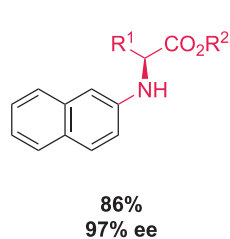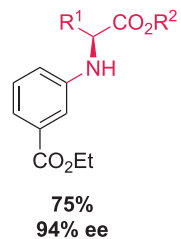


**Alcohol and Hydroxide Coupling<sup>1</sup>**

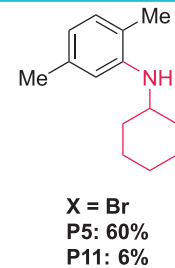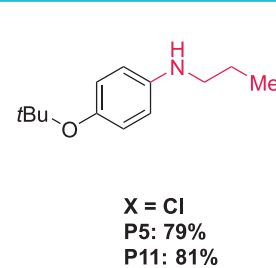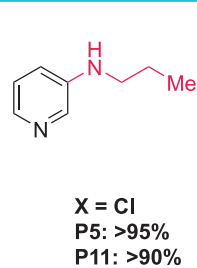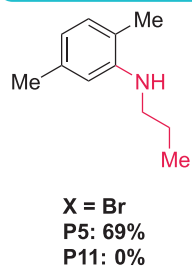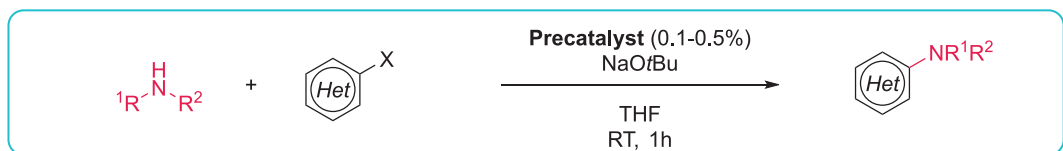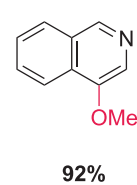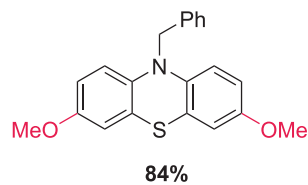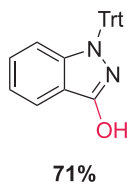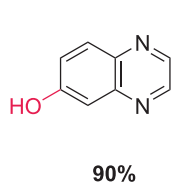


**Buchwald Hartwig Aminations with Primary Aliphatic Amines<sup>2</sup>**

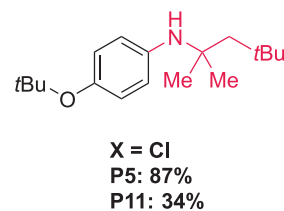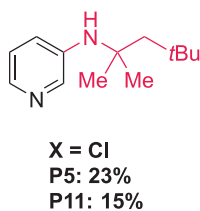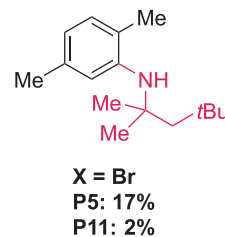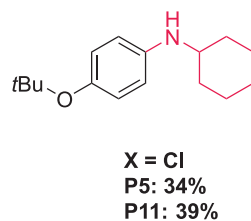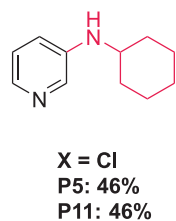


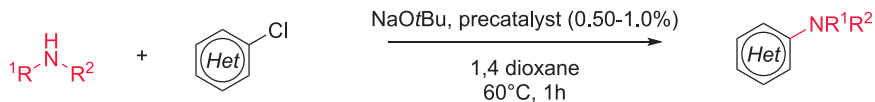


**Buchwald Hartwig  
Amination with  
Alkyl Amines and  
N Heterocycles<sup>3</sup>**

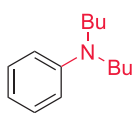


P5: 1%  
P12: 10%  
P14: 55%  
P15: 71%

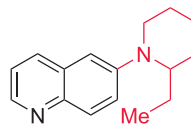


P5: 24%  
P12: 0%  
P14: 57%  
P15: 100%

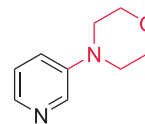


P5: 21%  
P12: 7%  
P14: 95%  
P15: 73%

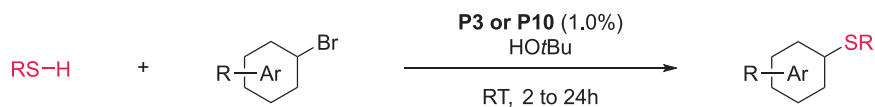


**Aliphatic  
Thiol Coupling  
of Hetero(aryl)  
Bromides<sup>4</sup>**

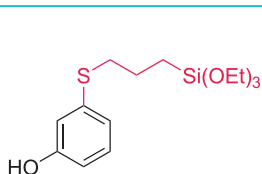


P3: 98%

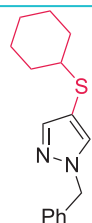


P3: 96%

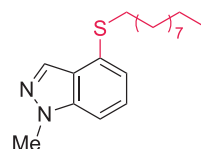

P10: 97%

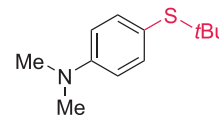


P10: 99%

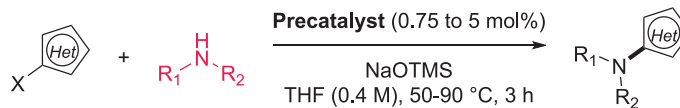


X = Br  
P4: 58%  
P5: 91%  
P6: 16%  
P11: 53%  
P12: 0%


**Buchwald-Hartwig  
Amination of  
Base-Sensitive Five-  
Membered Heteroaryl  
Halides and Aliphatic  
Amines<sup>5</sup>**

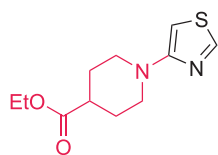


X = Br  
P5: 88%

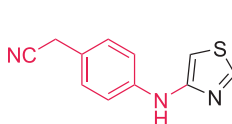


X = Cl  
P5: 90%

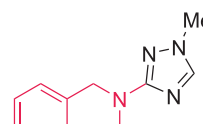


X = Br  
P5: 99%

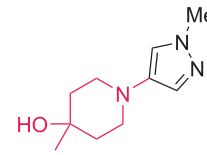


X = Br  
P5: 98%

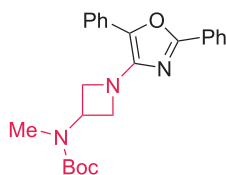


X = Br  
P5: 82%

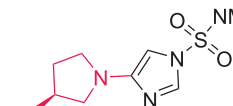

X = I  
P5: 94%

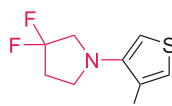

X = Br  
P5: 90%

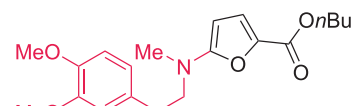

X = Br  
P5: 84%

**Table 2.** Buchwald G6 and Other OAC Precatalysts.

Precatalyst Reference ID	Product Number	Name	Structure
P1	915602	(AlPhos)Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	
P2	916455	(AlPhos)Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(OTf)	
P3	912883	( <i>t</i> BuBrettPhos)Pd( <i>p</i> -TMSCH <sub>2</sub> CH <sub>2</sub> CO <sub>2</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	
P4	937398	( <i>t</i> BuXPhos)Pd( <i>p</i> -MeC <sub>6</sub> H <sub>4</sub> )(Cl)	
P5	922900	(GPhos)Pd( <i>p</i> -TMSCH <sub>2</sub> CH <sub>2</sub> CO <sub>2</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	
P6	912646	( <i>t</i> BuBrettPhos)Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	

Precatalyst Reference ID	Product Number	Name	Structure
P7	915378	(AdBrettPhos)Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	
P8	925454	(AdCyBrettPhos)Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	
P9	931853	(SPhos)Pd( <i>p</i> -C <sub>4</sub> H <sub>4</sub> NO <sub>2</sub> CO <sub>2</sub> CH <sub>2</sub> CH <sub>2</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	
P10	936103	( <i>t</i> BuXPhos)Pd( <i>p</i> -C <sub>6</sub> H <sub>5</sub> OC(O)NHC <sub>6</sub> H <sub>4</sub> )(Br)	
P11	938734	(EPhos)Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	
P12	938742	(RuPhos)Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	
P13	944319	( <i>t</i> BuSPhos)Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	

Precatalyst Reference ID	Product Number	Name	Structure
P14	945625*	(3,5- <i>t</i> Bu-CyFPhos)Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	
P15	946230*	CyFPhos Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	
P16	937452*	(( <i>t</i> Bu)PhCPhos)Pd( <i>p</i> -CF <sub>3</sub> C <sub>6</sub> H <sub>4</sub> )(Br)	

\*Coming Soon

## Reference

- Ingoglia BT, Buchwald SL. 2017. Oxidative Addition Complexes as Precatalysts for Cross-Coupling Reactions Requiring Extremely Bulky Biarylphosphine Ligands. *Org. Lett.* 19(11):2853-2856. <https://doi.org/10.1021/acs.orglett.7b01082>
- McCann SD, Reichert EC, Arrechea PL, Buchwald SL. 2020. Development of an Aryl Amination Catalyst with Broad Scope Guided by Consideration of Catalyst Stability. *J. Am. Chem. Soc.* 142(35):15027-15037. <https://doi.org/10.1021/jacs.0c06139>
- Feng K, Reichert E, Raguram J, Howard JR, Peters E, Liu C, Sigman MS, Buchwald SL. 2024. Development of a Deactivation-Resistant Dialkylbiarylphosphine Ligand for Pd-Catalyzed Arylation of Secondary Amines. *J. Am. Chem. Soc.* 146(39):26609-26615. <https://doi.org/10.1021/jacs.4c09667>
- Xu J, Liu RY, Yeung CS, Buchwald SL. 2019. Monophosphine Ligands Promote Pd-Catalyzed C-S Cross-Coupling Reactions at Room Temperature with Soluble Bases. *ACS Catal.* 9(7):6461-6466. <https://doi.org/10.1021/acscatal.9b01913>
- Reichert EC, Feng K, Sather AC, Buchwald SL. 2023. Pd-Catalyzed Amination of Base-Sensitive Five-Membered Heteroaryl Halides with Aliphatic Amines. *J. Am. Chem. Soc.* 145(6):3323-3329. <https://doi.org/10.1021/jacs.2c13520>

To view our complete offerings, visit [SigmaAldrich.com/Buchwald](https://www.sigmaaldrich.com/buchwald)



To place an order or receive technical assistance:  
[SigmaAldrich.com/support](https://www.sigmaaldrich.com/support)



For local contact information:  
[SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices)

Merck KGaA  
Frankfurter Strasse 250  
64293 Darmstadt, Germany

[SigmaAldrich.com](https://www.sigmaaldrich.com)

We have built a unique collection of life science brands with unrivalled experience in supporting your scientific advancements.

**Millipore**® **Sigma-Aldrich**® **Supelco**® **Milli-Q**® **SAFC**® **BioReliance**®

© 2026 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. Merck, the vibrant M, BioReliance, Millipore, Milli-Q, SAFC, Sigma-Aldrich, and Supelco are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

Ver. 1.0 67656 03/2026

