(English version see page 2)
(Version français voir page 3)
(Versión en español vea página 4)

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Readme File for Firmware Update of Nova 30/60 **(firmware version 3.15 or higher)**

General information:

The program "Firmware Update Nova 30/60 " enables you to **update the firmware of the Nova 30/60** to the latest version with the aid of a PC. To do so, you need a free port on your PC and a suitable interface cable (see detailed operating manual on CD-ROM or Internet).

Program installation:

With the installation program "Nova_3060_Vx_yy.exe" you install the firmware update program on your PC.

For Microsoft Windows 7 or higher, administrator rights are required.

Program start:

In the Windows start menu, execute the program "Update Nova 3060" in the "Merck/Update Nova 3060" directory.

Via the Language menu, you can change the language (German or English).

For Microsoft Windows 7 or higher, administrator rights are required.

Firmware update:

Using a suitable cable (see detailed operating manual on CD-ROM or Internet), connect the Nova 30/60 to be programmed to an interface of the PC.

Make sure the instrument to be programmed is switched on.

Click the OK button to start the update process.

The program automatically recognizes the instrument connected.

Proceed by following the instructions of the program.

The programming process takes about 5 minutes. After successful programming, a final message appears. The programming process is finished now. Disconnect the instrument from the PC. The instrument is ready for operation. After switching Off/On you can check whether the new software version was taken over (menu setup/meter setup/system info).

Note:

Instrument functions can change with a firmware update. The current analysis specifications are available on the Internet under www.analytical-test-kits.com.

The current operating manual is available on the Internet under www.merckmillipore.com/photometry.

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