

# Product Information

## EX-CELL™ CD CHO Serum-Free Medium for CHO Cells, Chemically Defined

without L-glutamine, without hypoxanthine, without thymidine

CATALOG NO. 14360C

### Description

EX-CELL™ CD CHO is a chemically defined, animal-component free, serum-free liquid medium developed for the long-term growth of Chinese Hamster Ovary (CHO) cells and expression of antibodies or protein products in suspension culture. CHO suspension cultures can be subcultured directly into EX-CELL™ CD CHO from serum-supplemented or serum-free medium with little or no adaptation. EX-CELL™ CD CHO (Catalog No. 14360C) is formulated without hypoxanthine, thymidine and L-glutamine, making it an appropriate medium for selection systems such as DHFR<sup>+</sup> and Glutamine Synthetase (GS System™). For applications that do not require the selective pressure of a hypoxanthine/thymidine (HT)-deficient medium, we recommend the use of EX-CELL™ CD CHO with hypoxanthine and thymidine (Catalog No. 14361C).

### Formulation

The formula for EX-CELL™ CD CHO is proprietary to SAFC Biosciences. For additional information, please call our Technical Services department.

### Precautions

Use aseptic technique when handling or supplementing this medium. This product is for research or for further manufacturing use. THIS PRODUCT IS NOT INTENDED FOR HUMAN OR THERAPEUTIC USE.

### Storage

Store liquid medium at 2 to 8 C, protected from light. Do not use after the expiration date.

### Indications of Deterioration

Medium should be clear and free of particulates and flocculent material. Do not use if liquid medium is cloudy or contains precipitate. Other evidence of deterioration may include color change, pH shift or degradation of physical or performance characteristics.

### Preparation Instructions

EX-CELL™ CD CHO is formulated without L-glutamine. For applications requiring the use of L-glutamine, supplement with 8 mM L-glutamine by adding 40 mL/L of a 200 mM solution (Catalog No. 59202C). SAFC Biosciences recommends L-glutamine supplementation of the working volume only. Supplements, such as antibiotics, can be added to the sterilized medium using aseptic technique. Storage conditions and shelf life of the supplemented product may be affected by the nature of the supplement.

### Methods for Use

#### Adaptation

CHO cells that have been grown in attachment or suspension cultures in a conventional serum-supplemented medium can be readily grown in EX-CELL™ CD CHO with little or no adaptation. Adaptation to EX-CELL™ CD CHO requires healthy, viable cultures in mid-logarithmic growth phase. During adaptation, growth rates can be somewhat slower than normal expected results.

#### United States

SAFC Biosciences, Inc.  
13804 W. 107th Street  
Lenexa, Kansas 66215  
USA  
Phone +1 913-469-5580  
Toll free-USA 1 800-255-6032  
Fax +1 913-469-5584  
E-mail info-na@sial.com

#### Europe

SAFC Biosciences Ltd.  
Smeaton Road, West Portway  
Andover, Hampshire SP10 3LF  
UNITED KINGDOM  
Phone +44 (0)1264-333311  
Fax +44 (0)1264-332412  
E-mail info-eu@sial.com

#### Asia Pacific

SAFC Biosciences Pty. Ltd.  
18-20 Export Drive  
Brooklyn, Victoria 3025  
AUSTRALIA  
Phone +61 (0)3-9362-4500  
Toll free-AUS 1 800-200-404  
Fax +61 (0)3-9315-1656  
E-mail info-ap@sial.com

#### *Adaptation from attachment cultures*

1. Subculture the cells from serum-supplemented medium to EX-CELL™ CD CHO supplemented with 8 mM L-glutamine using standard trypsinization techniques when cultures reach 100% confluence.
2. Inactivate the trypsin with medium containing 5% gamma irradiated Fetal Bovine Serum (FBS) (Catalog No. 12107C). Using low-speed centrifugation, pellet the cell suspension at 200 *g* for 5 minutes and carefully decant the supernatant without disturbing the cell pellet.
3. Resuspend the cells in EX-CELL™ CD CHO medium supplemented with 8 mM L-glutamine at a density of  $5 \times 10^5$  cells/mL in 25 cm<sup>2</sup> or 75 cm<sup>2</sup> flasks.
4. Incubate the cells at 37 C in a humidified incubator with 8 - 10% CO<sub>2</sub>.
5. Continue to subculture the cells in EX-CELL™ CD CHO every 3 - 4 days for 3 - 4 passages using a seeding density of  $5 \times 10^5$  cells/mL in 25 cm<sup>2</sup> or 75 cm<sup>2</sup> flasks. SAFC Biosciences recommends centrifugation of the cell suspension prior to passaging.
6. After 3 - 4 passages, centrifuge and resuspend the cells in EX-CELL™ CD CHO medium supplemented with 8 mM L-glutamine at a density of  $3\text{-}5 \times 10^5$  cells/mL in 125 mL shaker flasks.
7. Continue to subculture cells in EX-CELL™ CD CHO every 3 - 4 days as described in step No. 6.
8. Allow the cells to adapt to EX-CELL™ CD CHO for an additional 4 - 6 passages. Cells are considered fully adapted to EX-CELL™ CD CHO when growth rates return to normal and viabilities are above 95%.
9. Cells can now be scaled up as necessary using standard procedures and densities.

#### *Adaptation from suspension cultures*

1. Resuspend the cells in EX-CELL™ CD CHO medium supplemented with 8 mM L-glutamine at a density of  $5 \times 10^5$  cells/mL in 125 mL shaker flasks.
2. Incubate the cells at 37 C in a humidified incubator with 8 - 10% CO<sub>2</sub>.
3. Continue to subculture the cells in EX-CELL™ CD CHO every 3 - 4 days for 3 - 4 passages using a seeding density of  $5 \times 10^5$  cells/mL in 125 mL flasks. SAFC Biosciences recommends centrifugation of the cell suspension prior to passaging.
4. After 3 - 4 passages, centrifuge and resuspend the cells in EX-CELL™ CD CHO medium supplemented with 8 mM L-glutamine at a density of  $3\text{-}5 \times 10^5$  cells/mL in 125 mL shaker flasks.

5. Continue to subculture cells in EX-CELL™ CD CHO every 3 - 4 days as described in step No. 4.
6. Allow the cells to adapt to EX-CELL™ CD CHO for an additional 4 - 6 passages. Cells are considered fully adapted to EX-CELL™ CD CHO when growth rates return to normal and viabilities are above 95%.
7. Cells can now be scaled up as necessary using standard procedures and densities.

#### **Culture Techniques**

CHO cells are normally grown at  $37 \pm 1$  C and 5 - 10% CO<sub>2</sub>. Allow the medium to warm to room temperature prior to use (protect from light). Once fully adapted, the cells should be subcultured at a seeding density of at least  $3 \times 10^5$  cells/mL in shaker flasks. Seed 30 mL cell cultures in 125 mL shaker flasks and 60 mL cultures in 250 mL shaker flasks. Shaker speed should be approximately 125 rpm  $\pm$  5 rpm. Optimal seeding densities and L-glutamine concentrations should be determined by the researcher for each application and cell type.

When passing the cells, medium carry over should not exceed 25% of the final volume. If carry over exceeds 25%, centrifugation is recommended. For best results, centrifugation is recommended regardless of medium carryover. Cells propagated in serum-free media are extremely fragile. For successful results, care must be taken when subculturing cells. Standard techniques of centrifugation must be modified to include low-speed centrifugation to prevent damage to cells that have been propagated in serum-free medium.

During adaptation, normal trypsin concentrations may be used, but incubations should be carried out at 4 C, and exposure time should be minimal. SAFC Biosciences recommends the use of a soybean trypsin inhibitor (0.1%), or sedimentation by centrifugation to remove the trypsin. Soybean trypsin inhibitor should be used with caution, as it is toxic to some cell types. Cells may also be dislodged with NO-ZYME™ (Catalog No. 59226C), a non-enzymatic dissociating agent.

## Cryopreservation

### *Freezing:*

Cells can be frozen in EX-CELL™ CD CHO without the reintroduction of serum. However, it is necessary to handle the cells gently and freeze the cells under carefully controlled conditions.

1. Choose cultures in logarithmic growth with viabilities above 90%.
2. Prepare a freezing medium consisting of 45% cold EX-CELL™ CD CHO medium, 45% conditioned medium and 10% dimethyl sulfoxide (DMSO). Alternatively, cells can be frozen in 90% fresh medium and 10% DMSO.
3. Centrifuge the cells at 200 *g* for 5 minutes. Remove the supernatant.
4. Resuspend the cells in the freezing medium at  $1 \times 10^7$  cells/mL.
5. Rapidly transfer 1 - 2 mL of this suspension to sterile cryovials.
6. Place the vials at -20 C for 3 - 4 hours, then transfer to -70 C for 16 - 24 hours.
7. For long-term storage, transfer the vials to liquid nitrogen vapor.

### *Thawing:*

1. Rapidly thaw a vial of frozen cells in a 37 C water bath.
2. Transfer the cells aseptically to a centrifuge tube containing 5 - 10 mL of cold EX-CELL™ CD CHO medium.
3. Using low-speed centrifugation, pellet the cell suspension at 200 *g* for 5 minutes and carefully decant the supernatant without disturbing the cell pellet.
4. Resuspend the cells in 5 mL of EX-CELL™ CD CHO.
5. Count the cells for viability and transfer to a sterile shaker flask at a seeding density of  $4 \times 10^5$  cells/mL.

## Characteristics

### **Appearance**

Clear yellow solution

### **Endotoxin**

Refer to Certificate of Analysis

### **Osmolality (as supplied)**

310 - 350 mOsm/kg H<sub>2</sub>O

### **pH (as supplied)**

7.0 - 7.4

### **Sterility**

No microbial growth detected

#### Warranty, Limitation of Remedies

SAFC Biosciences warrants to the purchaser for a period of one year from date of delivery that this product conforms to its specifications. Other terms and conditions of this warranty are contained in SAFC Biosciences' written warranty, a copy of which is available upon request. ALL OTHER WARRANTIES, EXPRESSED OR IMPLIED, INCLUDING THE IMPLIED WARRANTY OF MERCHANTABILITY AND FITNESS FOR A PARTICULAR PURPOSE, ARE EXCLUDED. In no case will SAFC Biosciences be liable for any special, incidental, or consequential damages arising out of this product or the use of this product by the customer or any third party based upon breach of warranty, breach of contract, negligence, strict tort, or any other legal theory. SAFC Biosciences expressly disclaims any warranty against claims by any third party by way of infringement or the like. THIS PRODUCT IS INTENDED FOR PURPOSES DESCRIBED ONLY AND IS NOT INTENDED FOR ANY HUMAN OR THERAPEUTIC USE.

Additional Terms and Conditions are contained in the product Catalog, a copy of which is available upon request.

EX-CELL™ and NO-ZYME™ are trademarks of SAFC Biosciences, Inc.  
The GS System™ is a trademark of Lonza Biologics.

© 2006 SAFC Biosciences, Inc.

Issued September 2006 P14360  
0204 0805 0905 0306

#### United States

SAFC Biosciences, Inc.  
13804 W. 107th Street  
Lenexa, Kansas 66215  
USA  
Phone +1 913-469-5580  
Toll free-USA 1 800-255-6032  
Fax +1 913-469-5584  
E-mail info-na@sial.com

#### Europe

SAFC Biosciences Ltd.  
Smeaton Road, West Portway  
Andover, Hampshire SP10 3LF  
UNITED KINGDOM  
Phone +44 (0)1264-333311  
Fax +44 (0)1264-332412  
E-mail info-eu@sial.com

#### Asia Pacific

SAFC Biosciences Pty. Ltd.  
18-20 Export Drive  
Brooklyn, Victoria 3025  
AUSTRALIA  
Phone +61 (0)3-9362-4500  
Toll free-AUS 1 800-200-404  
Fax +61 (0)3-9315-1656  
E-mail info-ap@sial.com

**www.safcbiosciences.com**