

17774 Potassium tellurite Solution

A filter sterilized solution of potassium tellurite, recommended for the selective isolation of staphylococci and corneobacteria.

Composition

(10 ml per vial)

sterile 1% Potassium tellurite Solution in water

Directions

Warm up potassium tellurite (1%) to 45-50°C and use in an appropriate agar or broth as required.

Supplement for: Vogel-Johnson Agar 70195 (10 ml per 500 ml), Baird Parker Agar, Base 11705 (10.5 ml per litre), Giolitti-Cantoni-Broth (0.1-0.01%), Mitis Salivarius Agar 01337 (1 ml per litre).

Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

