

## Technical Data Sheet

# KF Streptococcus Agar (base)

Ordering number: 1.10707.0500

For the detection and enumeration of enterococci (fecal streptococci) in water, foodstuffs and other materials according to KENNER, CLARK and KABLER (1960, 1961).

KF (Kenner Fecal) Streptococcus agar complies with the recommendations given by APHA for the examination of water (1998) and foodstuffs (1992).

### Mode of Action

Maltose and lactose are metabolized by most enterococci with the production of acid and thus promote the growth of these bacteria; undesired microorganisms are largely suppressed by sodium azide. Acid formation is detected by bromocresol purple which changes its colour to yellow. Enterococci reduce TTC to give a red formazan and thus appear as red colonies.

### Typical Composition (g/L)

| KF Streptococcus Agar Base |       |
|----------------------------|-------|
| Proteose peptone           | 10.0  |
| Yeast extract              | 10.0  |
| Sodium chloride            | 5.0   |
| Sodium glycerophosphate    | 10.0  |
| Maltose                    | 20.0  |
| Lactose                    | 1.0   |
| Sodium azide               | 0.4   |
| Bromocresol purple         | 0.015 |
| Agar-agar**                | 15.0  |

\*\*Agar-agar is equivalent to other different terms of agar.

### Also to be added:

2,3,5-triphenyltetrazolium chloride 0.1

## Preparation

Suspend 71.5 g in 1 litre of demineralised water. Boil up with frequent agitation. Boil for 5 minutes (or autoclave 10 min at 121°C, if total selectivity is required).

**Do not overheat.**

Cool to approx. 50 °C, add 10 ml of a 1 % TTC solution (2,3,5- triphenyltetrazolium chloride), mix, pour plates.

pH: 7.2 ± 0.2 at 25 °C.

The plates are clear and purple.

## Experimental Procedure and Evaluation

The membrane filtration method should be used for detection and enumeration if only small numbers of enterococci are suspected to be present; the pour plate method should be employed for larger numbers. The inoculated membrane filters are placed on the agar surface.

Incubation: 48 hours at 35 °C aerobically.

Tropical marine water samples should be incubated anaerobically, due to high incidence of false-positive presumptive counts for enterococci.

The red or pink colonies should be counted, the bacterial count can then be calculated.

| Appearance of Colonies                                            | Microorganisms                                                                                                                                                  |
|-------------------------------------------------------------------|-----------------------------------------------------------------------------------------------------------------------------------------------------------------|
| Abundant growth, red colonies, mostly surrounded by a yellow zone | Enterococci (E. faecalis, E. faecalis var. liquefaciens, E. faecalis var. zymogenes), Streptococcus mitis, Str. bovinus, E. equinus, Str. salivarius and others |
| Usually scanty growth and no colour change                        | Lactobacillus plantarum, Pediococcus cerevisiae and others                                                                                                      |

## Storage

Store at +15 °C to +25 °C, dry and tightly closed. Do not use clumped or discolored medium. Protect from UV light (including sun light). For *in vitro* use only.



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
We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

Merck Millipore, the M logo, are registered trademarks of Merck KGaA, Darmstadt, Germany. ATCC is a registered trademark of ATCC, Manassas, VA, USA. Lit. No. TN1822EN00

## Quality Control

| Control strains                               | Growth, 48h, 35°C, aerobic | Red colonies | Colour change to yellow |
|-----------------------------------------------|----------------------------|--------------|-------------------------|
| Enterococcus faecalis ATCC 11700              | Good to very good          | +            | +                       |
| Enterococcus faecalis ATCC 19433 (WDCM 00009) | Good to very good          | +            | +                       |
| Enterococcus hirae ATCC 8043 (WDCM 00089)     | Good to very good          | +(poor)      | +                       |
| Streptococcus pyogenes ATCC 12344             | None to fair               | -            | -                       |
| Streptococcus agalactiae ATCC 13813           | None to fair               | -            | -                       |
| Lactobacillus plantarum ATCC 8014             | None to fair               | -            | -                       |
| Escherichia coli ATCC 25922 (WDCM 00013)      | None                       |              |                         |
| Enterobacter cloacae ATCC 13047               | None                       |              |                         |
| Pseudomonas aeruginosa ATCC 27853             | none                       |              |                         |

Please refer to the actual batch related Certificate of Analysis.



Enterococcus faecalis ATCC 11700



Streptococcus agalactiae ATCC 13813



We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

Merck Millipore, the M logo, are registered trademarks of Merck KGaA, Darmstadt, Germany. ATCC is a registered trademark of ATCC, Manassas, VA, USA. Lit. No. TN1822EN00

## Literature

American Public Health Association: Compendium of methods for the microbiological examination of foods. – 3 rd. ed., 1992.

American Public Health Association: American Water Works Association and Water Pollution Control Federation: Standard Methods for the Examination of Water and Wastewater 20 th ed., Washington, 1998.

KENNER, B.A., CLARK, H.F., a. KABLER F.W.: Faecal streptococci. II. Quantification of streptococci in faeces. - **Am. J. Publ. Health.**, **50**; 1553-1559 (1960).

KENNER, B.A., CLARK, H.F., a. KABLER F.W.: Faecal streptococci. I. Cultivation and enumeration of streptococci in surface waters. - **Appl. Microbiol.**, **9**; 15-20 (1961)

## Ordering Information

| Product                             | Cat. No.      | Pack size |
|-------------------------------------|---------------|-----------|
| KF Streptococcus Agar Base          | 1.10707 .0500 | 500 g     |
| 2,3,5-Triphenyltetrazolium chloride | 1.08380.0010  | 10 g      |

Merck KGaA, 64271 Darmstadt, Germany  
Fax: +49 (0) 61 51 / 72-60 80  
mibio@merckgroup.com  
[www.merckmillipore.com/biomonitoring](http://www.merckmillipore.com/biomonitoring)

Find contact information for your country at:  
[www.merckmillipore.com/offices](http://www.merckmillipore.com/offices)  
For Technical Service, please visit:  
[www.merckmillipore.com/techservice](http://www.merckmillipore.com/techservice)



We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

Merck Millipore, the M logo, are registered trademarks of Merck KGaA, Darmstadt, Germany. ATCC is a registered trademark of ATCC, Manassas, VA, USA. Lit. No. TN1822EN00