



SIGMA-ALDRICH

3050 Spruce Street
Saint Louis, Missouri 63103 USA
Telephone 800-325-5832 • (314) 771-5765
Fax (314) 286-7828
email: techserv@sial.com
sigma-aldrich.com

ProductInformation

SIGMA QUALITY CONTROL TEST PROCEDURE

Enzymatic Assay of TRYPSIN¹ (EC 3.4.21.4)

PRINCIPLE:



Abbreviation used:

BAEE = N α -Benzoyl-L-Arginine Ethyl Ester

CONDITIONS: T = 25EC, pH = 7.6, A_{253nm}, Light path = 1 cm

METHOD: Continuous Spectrophotometric Rate Determination

REAGENTS:

- A. 67 mM Sodium Phosphate Buffer, pH 7.6 at 25EC
(Prepare 100 ml in deionized water using Sodium Phosphate, Monobasic, Anhydrous, Sigma Prod. No. S-0751. Adjust to pH 7.6 at 25EC with 1 M NaOH.)
- B. 0.25 mM N α -Benzoyl-L-Arginine Ethyl Ester Solution (BAEE)
(Prepare 50 ml in Reagent A using N α -Benzoyl-L-Arginine Ethyl Ester, Hydrochloride, Sigma Prod. No. B-4500.)
- C. 1 mM Hydrochloric Acid Solution (HCl)
(Prepare 50 ml in deionized water using concentrated Hydrochloric Acid, Sigma Prod. No. H-7020.)
- D. Trypsin Enzyme Solution
(Immediately before use, prepare a solution containing 500 BAEE units/ml of Trypsin in cold Reagent C.)

Enzymatic Assay of TRYPSIN¹ (EC 3.4.21.4)

PROCEDURE:

Pipette (in milliliters) the following reagents into suitable quartz cuvettes:

	<u>Test</u>	<u>Blank</u>
Reagent B (BAEE)	3.00	3.00

Equilibrate to 25EC. Monitor the A_{253nm} until constant, using a suitably thermostatted spectrophotometer. Then add:

Reagent C (HCl)	-----	0.20
Reagent D (Enzyme Solution)	0.20	-----

Immediately mix by inversion and record the increase in A_{253nm} for approximately 5 minutes. Obtain the $\Delta A_{253nm}/\text{minute}$ using the maximum linear rate for both the Test and Blank.

CALCULATIONS:

$$\text{BAEE Units/ml enzyme} = \frac{(\Delta A_{253nm}/\text{min Test} - \Delta A_{253nm}/\text{min Blank})(df)}{(0.001) (0.20)}$$

df = Dilution factor

0.001 = The change in A_{253nm}/minute per unit of Trypsin at pH 7.6 at 25EC in a 3.2 ml reaction mix

0.20 = Volume (in milliliters) of enzyme used

$$\text{Units/mg solid} = \frac{\text{units/ml enzyme}}{\text{mg solid/ml enzyme}}$$

$$\text{Units/mg protein} = \frac{\text{units/ml enzyme}}{\text{mg protein/ml enzyme}}$$

UNIT DEFINITION:

One BAEE unit will produce a ΔA_{253nm} of 0.001 per minute with BAEE as substrate at pH 7.6 at 25EC in a reaction volume of 3.2 ml.

FINAL ASSAY CONCENTRATION:

In a 3.2 ml reaction mix, the final concentrations are 63 mM sodium phosphate, 0.23 mM N α -benzoyl-L-arginine ethyl ester, 0.06 mM hydrochloric acid, and 100 units trypsin.

Enzymatic Assay of TRYPSIN¹ **(EC 3.4.21.4)**

REFERENCE:

Bergmeyer, H.U., Gawehn, K., and Grassl, M. (1974) in *Methods of Enzymatic Analysis* (Bergmeyer, H.U., ed) Volume I, 2nd ed., 515-516, Academic Press, Inc., New York, NY

NOTES:

1. This assay procedure is not to be used to assay Sigma Prod. Nos. T-1763, T-4019, T-8386, T-8899, and T-9906.
2. This assay is based on the cited reference.
3. Where Sigma Product or Stock numbers are specified, equivalent reagents may be substituted.

Sigma warrants that the above procedure information is currently utilized at Sigma and that Sigma products conform to the information in Sigma publications. Purchaser must determine the suitability of the information and products for its particular use. Upon purchase of Sigma products, see reverse side of invoice or packing slip for additional terms and conditions of sale.