

1.09874.0010

Microbiology

MSRV Selective Supplement

Contents: 10 vials of lyophilisate for the preparation of 5.0 l
MSRV Medium

Composition (per vial)

Novobiocin 10 mg

Preparation of MSRV Medium

1. Suspend 15.8 g MSRV Medium (Base) in 500 ml of demin. water by heating in a boiling water bath or in a flowing steam until the medium is completely dissolved.
Do not autoclave!
2. Let the medium cool to 45–50 °C.
3. Dissolve the supplement by adding 1 ml sterile distilled water and add the solution to the medium. Mix gently and pour to plates.

Ordering information

Cat. No.	Product	Package size
1.09878.0500	MSRV Medium (Base)	500 g

Status: 2009-10-19

Merck KGaA
64271 Darmstadt, Germany
Tel. + 49 (0)6151 72-2440
www.merck-chemicals.com

