


Anti-mTOR (T 2949)

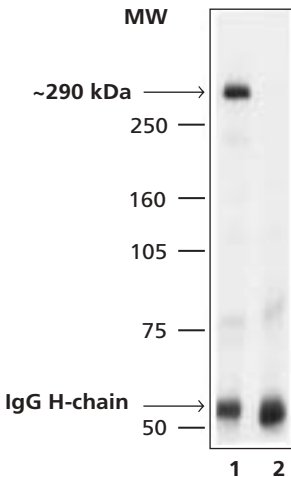

Immunoprecipitation

mTOR was immunoprecipitated from whole extract of human HEK 293T cells using Rabbit Anti-mTOR (Product No. T 2949). The IP samples were separated on SDS-PAGE and blotted with a control Rabbit anti-rat mTOR antibody. The antibody was developed with Goat Anti-Rabbit IgG, Peroxidase Conjugate (Product No. A 0545) and chemiluminescent substrate.

Lane 1: 1.1 μ g of Anti-mTOR immunoprecipitates mTOR from a RIPA extract of 300 μ g of 293T cells.

Lane 2: negative control (pre-immune serum was used instead of the Anti-mTOR antibody)