


## Product Information

### Sigma-Aldrich 06929 Timestrip® 12 Months, Key Chain




The Timestrip® 12 Months key chain indicator monitors time elapsed over a period of twelve months. It is designed to be used at room temperature conditions. It is inert until activated and can be stored at room temperature until needed. Housed in a handy key chain, it is self-adhesive, single use and irreversible. It has a pre-activation shelf life of three years from shipping.


#### Instruction for use

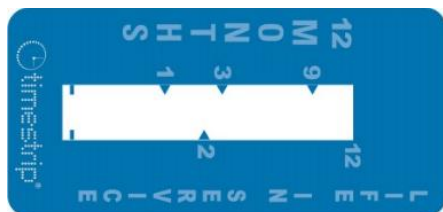

**To Activate:** Remove the small stick on the key chain and push the black lever down into the clear plastic holder. This will press the button on the back of the indicator and release the red dye. A red line will appear to show that the indicator has been activated.

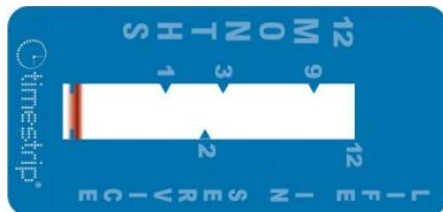

**Activated:** Attach the indicator in the key chain to the product or device to be monitored.

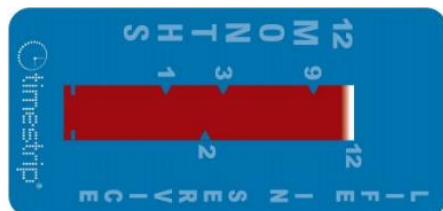

**Review:** The red dye will move through the viewing window and past the time markers. When the red dye reaches the end of the viewing window, twelve months have passed.

For full instructions, testing protocols, specifications and other information on this product, please visit [www.sigmaaldrich.com/timestrip](http://www.sigmaaldrich.com/timestrip)