




Product Information

SIGMA QUALITY CONTROL TEST PROCEDURE

Enzymatic Assay of NITRATE REDUCTASE (NAD[P]H) (EC 1.6.6.2) Sigma Prod. No. N-7265

PRINCIPLE:

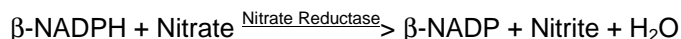

Abbreviations used:

β -NADPH = β -Nicotinamide Adenine Dinucleotide Phosphate, Reduced Form

β -NADP = β -Nicotinamide Adenine Dinucleotide Phosphate, Oxidized Form

CONDITIONS: T = 25°C, pH = 7.5, $A_{340\text{nm}}$, Light path = 1 cm

METHOD: Continuous Spectrophotometric Rate Determination

REAGENTS:

- A. 100 mM Potassium Phosphate Solution
(Prepare 100 ml in deionized water using Potassium Phosphate, Monobasic, Anhydrous, Sigma Prod. No. P-5379.)
- B. 100 mM Potassium Phosphate Buffer, pH 7.5 at 25°C
(Prepare 100 ml in deionized water using Potassium Phosphate, Dibasic, Trihydrate, Sigma Prod. No. P-5504. Adjust to pH 7.5 at 25°C with Reagent A.)
- C. 0.1 mM Flavin Adenine Dinucleotide Solution (FAD)
(Prepare 100 ml in deionized water using Flavin Adenine Dinucleotide, Disodium Salt, Sigma Prod. No. F-6625. **PREPARE FRESH and PROTECT FROM LIGHT.**)
- D. 12 mM β -Nicotinamide Adenine Dinucleotide Phosphate, Reduced Form Solution (β -NADPH)
(Dissolve the contents of one 10 mg vial of β -Nicotinamide Adenine Dinucleotide Phosphate, Reduced Form, Tetrasodium Salt, Sigma Stock No. 201-210, in the appropriate volume of deionized water. **PREPARE FRESH.**)
- E. 100 mM Potassium Nitrate Solution (KNO_3)
(Prepare 1 ml in deionized water using Potassium Nitrate, Sigma Prod. No. P-8394.)

**Enzymatic Assay of NITRATE REDUCTASE (NAD[P]H)
(EC 1.6.6.2)
Sigma Prod. No. N-7265**

REAGENTS: (continued)

- F. Nitrate Reductase Enzyme Solution
(Immediately before use, prepare a solution containing 0.13 - 0.25 unit/ml of Nitrate Reductase in cold Reagent B. Store on ice; use within 5 minutes.)

PROCEDURE:

Pipette (in milliliters) the following reagents into suitable cuvettes:

	<u>Test</u>	<u>Blank</u>
Deionized Water	0.80	1.10
Reagent B (Buffer)	1.60	1.60
Reagent C (FAD)	0.15	0.15
Reagent D (β -NADPH)	0.05	0.05
Reagent E (KNO_3)	0.30	-----

Mix by inversion and equilibrate to 25°C. Monitor the $A_{340\text{nm}}$ until constant, using a suitably thermostatted spectrophotometer. Then add:

Reagent F (Enzyme Solution)	0.10	0.10
-----------------------------	------	------

Immediately mix by inversion and record the decrease in $A_{340\text{nm}}$ for approximately 5 minutes. Obtain the $\Delta A_{340\text{nm}}$ /minute using the maximum linear rate for both the Test and Blank.

CALCULATIONS:

$$\text{Units/ml enzyme} = \frac{(\Delta A_{340\text{nm}}/\text{min Test} - \Delta A_{340\text{nm}}/\text{min Blank})(3)(\text{df})}{(6.22)(0.1)}$$

3 = Total volume (in milliliters) of assay

df = Dilution factor

6.22 = Millimolar extinction coefficient of β -NADPH at 340 nm

0.1 = Volume (in milliliter) of enzyme used

$$\text{Units/mg solid} = \frac{\text{units/ml enzyme}}{\text{mg solid/ml enzyme}}$$

**Enzymatic Assay of NITRATE REDUCTASE (NAD[P]H)
(EC 1.6.6.2)
Sigma Prod. No. N-7265**

CALCULATIONS: (continued)

$$\text{Units/mg protein} = \frac{\text{units/ml enzyme}}{\text{mg protein/ml enzyme}}$$

UNIT DEFINITION:

One unit will reduce 1.0 μ mole of nitrate per minute in the presence of β -NADPH at pH 7.5 at 25°C.

FINAL ASSAY CONCENTRATION:

In a 3.00 ml reaction mix, the final concentrations are 57 mM potassium phosphate, 0.005 mM flavin adenine dinucleotide, 0.2 mM β -nicotinamide adenine dinucleotide phosphate, reduced form, 10 mM potassium nitrate and 0.013 - 0.025 unit nitrate reductase.

REFERENCE:

Gilliam, M.B., Sherman, M.P., Griscavage, J.M., and Ignarro, L.J. (1993) *Analytical Biochemistry* **212**, 359-365

NOTES:

1. This assay is based on the cited reference.
2. Where Sigma Product or Stock numbers are specified, equivalent reagents may be substituted.

Sigma warrants that the above procedure information is currently utilized at Sigma and that Sigma products conform to the information in Sigma publications. Purchaser must determine the suitability of the information and products for its particular use. Upon purchase of Sigma products, see reverse side of invoice or packing slip for additional terms and conditions of sale.