

## Product Information

# Adenosine 5'-monophosphate Agarose

Lyophilized powder

**A1271**

## Product Description

Adenosine 5'-monophosphate-Agarose (5'-AMP-agarose) is a conjugate of 5'-AMP to cross-linked 4% beaded agarose (activated by cyanogen bromide), via the C-8 atom of 5'-AMP. 5'-AMP-agarose is used in studies of NAD<sup>+</sup>-dependent dehydrogenases and ATP-dependent enzymes.

5'-AMP-agarose is useful in affinity chromatography for purifying various enzymes, such as:

- Aldehyde dehydrogenase<sup>1</sup>
- DNA polymerase  $\delta^2$
- DNA polymerase  $\epsilon^3$
- Lactate Dehydrogenase Isoenzyme-5<sup>4</sup>
- NADH oxidase<sup>5</sup>

## Storage/Stability

- The lyophilized 5'-AMP-agarose resin should be stored at -20 °C.
- Hydrated 5'-AMP-agarose resin can be stored refrigerated in neutral pH buffer that contains a bacteriostat, such as 0.02% sodium azide or thimerosal, or 20% ethanol.
- Do not autoclave or freeze the hydrated resin.
- The resin can be used several times without loss of effectiveness.

## Preparation Instructions

- The 5'-AMP-agarose resin may be hydrated by the addition of excess water, e.g. 100-150 mL per gram of resin, depending on the observed degree of swelling, for at least 30 minutes.
- The lactose stabilizer may be removed by washing the resin on a Buchner funnel with gentle vacuum, using 100-200 mL of water per gram of resin.
- Do not allow the resin to dry.
- The washed resin may then be resuspended in excess water or starting buffer, to pack the column bed.

## Procedure

Used 5'-AMP-agarose resin may be regenerated by washing the resin with 2-3 bed volumes of buffers in a cycle with alternating mildly basic and mildly acidic buffers, such as the following buffers:

- 0.1 M Trizma®-HCl plus 0.5 M NaCl, pH 8.5
- 0.1 M sodium acetate plus 0.5 M NaCl, pH 4.5

Three buffer wash cycles are suggested. The washed resin should then be re-equilibrated in 3-5 bed volumes of the desired binding buffer.

In situations where difficult-to-elute materials, such as lipids or denatured proteins, have become entrapped on the resin, a dilute detergent solution, such as 0.1% Triton™ X-100, may be used to wash the resin at 37 °C. The resin should then be immediately re-equilibrated with a minimum of 5 bed volumes of the desired binding buffer.

## References

1. Murphy, C.D. *et al.*, *Appl. Environ. Microbiol.*, **67(10)**, 4919-4921 (2001).
2. Cho, S.-W. *et al.*, *Mol. Cells*, **5(1)**, 20-24 (1995).
3. Lee, S.-K., and Fuchs, M.S., *FEBS Lett.*, **316(3)**, 261-263 (1993).
4. Pettit, S.M. *et al.*, *Clin. Chem.*, **27(1)**, 88-93 (1981).
5. Brown, D.M. *et al.*, *Eur. J. Biochem.*, **241(1)**, 155-161 (1996).

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

## Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice).

## Standard Warranty

The applicable warranty for the products listed in this publication may be found at [SigmaAldrich.com/terms](https://www.sigmaaldrich.com/terms).

## Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://www.sigmaaldrich.com/offices).

The life science business of Merck operates as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2021 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

A1271pis Rev 09/21 GCY,MAM

For research use only. Not for use in diagnostic procedures.

The Merck logo is displayed in a bold, red, sans-serif font.