Safety Sheet

mPAGE[®] Lux Curing Station

LUXGCS

Select language:

English	2	한국어 (Korean)	30
Български (Bulgarian)	4	Latviski (Latvian)	32
中文 (Chinese, Simplified) . .	6	Lietuvių (Lithuanian)	34
Hrvatski (Croatian)	8	Norsk (Norwegian)	36
Čeština (Czech)	10	Polski (Polish)	38
Dansk (Danish)	12	Português (Portuguese) . .	40
Engels (Dutch)	14	Română (Romanian)	42
Eesti keel (Estonian)	16	Srpski (Serbian)	44
Français (French)	18	Slovenčina (Slovak)	46
Deutsch (German)	20	Slovenščina (Slovenian) . .	48
Ελληνικά (Greek)	22	Español (Spanish)	50
Magyar (Hungarian)	24	Svenska (Swedish)	52
Italiano (Italian)	26	Türkçe (Turkish)	54
日本語 (Japanese)	28		

Intended Use

The mPAGE® Lux Curing Station emits ultraviolet (UV) light of a specified wavelength to initiate the polymerization of mPAGE® Lux Bis-Tris gel solutions to form a polyacrylamide gel. The gels produced are used for the separation of protein mixtures via electrophoresis, a process known as polyacrylamide gel electrophoresis (PAGE). The mPAGE® Lux Curing Station is intended only for use with mPAGE® Lux Bis-Tris Reagent Kits.

Specifications

Device

Rated Voltage	24 VDC		
Rated Power	65 W		
Dimensions	Height	20.32 cm	8 in.
	Width	26.04 cm	10.25 in.
	Depth	27.94 cm	11 in.
Weight	3.084 kg	6.8 lbs	

External Power Supply

Input Voltage	100-240 VAC (± 10%)
Input Current	1.5 A
Input Frequency	50/60 hz
Output Voltage	24 VDC
Output Current	2.7 A

Environmental Conditions

For indoor use only. This equipment is not rated for wet locations or any other outdoor parameter.

Storage

Temperature	15 °C to 30 °C
Humidity	20-80%, non-condensing

Operating

Temperature	15 °C to 30 °C
Humidity	20-80%, non-condensing
Atmospheric Pressure	Below 2000 M

For Indoor use only

Intended for Pollution Degree II
based on IEC61010-1

User Guide

User guides and other documents are available online. Go to [SigmaAldrich.com](https://www.sigmaaldrich.com), enter the catalogue number in the search bar, select your product, and scroll to the documents section.

The User Guide contains a list of consumables required to operate this device. The can be ordered online at [SigmaAldrich.com](https://www.sigmaaldrich.com).

Precautions

Please read these safety instructions carefully before using the device and/or any accessory supplied with this product. Keep a printed copy of the User Guide near the system for reference.

General Warnings

- This device complies with part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) This device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.
- Changes or modifications could void the user's authority to operate the equipment.
- During initial startup system will run self-test to insure device is functioning correctly.
- Ultraviolet (UV) light is emitted from this product. Minimize exposure to eyes and skin. Use appropriate shielding.
- Do not use with flammable or explosive liquids.
- Do not immerse device body in liquid.
- Do not expose device to vibrations. Vibrations may cause device malfunction or damage.
- Use only authorized accessories.
- If the device is broken or dropped, do not try to repair the device. Go to [SigmaAldrich.com/techservice](https://www.sigmaaldrich.com/techservice) for assistance.

Electrical Warnings

Failure to follow any of these warnings may result in injury, fire, and electric shock.

- Do not touch plug with wet hands.
- To avoid potential shock hazard, choose the correct plug configuration and ensure the power adapter is plugged securely into a properly grounded AC power outlet.
- Always ensure that the power supply input voltage matches the voltage available in your location.
- To avoid danger of electric shock, do not install the device in an area with a high humidity level, such as a greenhouse or an incubator. Refer to Environmental Conditions.
- To avoid damage to internal electronic components, do not spray device with sanitizing agents. Refer to Maintenance section in the User Guide for sanitizing information.
- Be sure that the mains power cord and plug are easily accessible in the case of a hazard in which the cord and plug need to be pulled in order to remove power from the product.

Maintenance

- Power OFF and unplug the instrument before beginning cleaning, to prevent risk of electrical shock.
- Do not spray any cleaner directly onto the instrument. **DO NOT ATTEMPT TO CLEAN ANY INTERNAL PARTS OF THE MACHINE.**
- Wipe clean with water. If needed, use mild detergent. Dry thoroughly before next use.
- Do not expose to acetone or ethyl alcohol. Do not use corrosive, flammable, or caustic cleaning agents.
- The instrument is NOT autoclave-able. Extreme heat will damage electronic components.
- If liquid enters the instrument, contact SigmaAldrich.com/techservice.

Technical Service

For assistance, go to SigmaAldrich.com/techservice. Unauthorized repair or service to the device may void the warranty and/or cause hazardous conditions.

Equipment Operation

See the user guide for protocols and use instructions.

Power ON

Locate the power button on the front panel of the device. Press the power button. The LED will light up above POWER, indicating power is ON.

Power OFF

Press and hold the power button for 5 seconds. The light above POWER will turn off, confirming power is off.

Cancel Cycle

Press "Start/Cancel" or open the door to cancel a cycle once initiated.

Emergency Disconnect

In an emergency, quickly unplug the device power cord from the mains power supply. Contact SigmaAldrich.com/techservice before restarting the device.

Power Button

Symbol Definitions

	Consult this Safety Sheet for important cautionary information
	Risk of injury by electrical shock
	UV light emitted from this product. Minimize exposure to eyes and skin by using appropriate shielding
	Catalogue Number
	Serial Number
	Lot Number
	Manufacturer

Declarations of Conformity are available on the product page at SigmaAldrich.com.

	This product has been evaluated by an OSHA certified Nationally Recognized Testing laboratory (NRTL) in the United States. Conforms to UL Std 61010-1 Certified to CSA Std 61010-1
	European Council Directives conformity marking.
	United Kingdom conformity marking

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Standard Warranty

The applicable warranty may be found at SigmaAldrich.com/terms.

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Made in USA

EMD Millipore Corporation,
Burlington, MA, U.S.A.,
an affiliate of Merck KGaA,
Darmstadt, Germany

MERCK

The life science business of Merck operates
as MilliporeSigma in the U.S. and Canada.

Merck, Millipore, mPAGE, and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

20723284w Rev 10/22