

Product Information

Sigma-Aldrich 07603 Timestrip® 3 Months



The Timestrip® 3 Months indicator monitors time elapsed over a period of three months. It is designed to be used at room temperature conditions. It is inert until activated and can be stored at room temperature until needed. It is self-adhesive, single use and irreversible. It has a pre-activation shelf life of three years from shipping.

Instruction for use



To Activate: Fully squeeze the button on top of the indicator. A red line will appear to show that the indicator has been activated.



Activated: Peel off the liner and stick the indicator to the product or device to be monitored.



Review: The red dye will move through the viewing window and past the time markers. When the red dye reaches the end of the viewing window, three months have passed.

For full instructions, testing protocols, specifications and other information on this product, please visit www.sigmaaldrich.com/timestrip