

Data Sheet

BioTracker™ Apo-15 Calcium-Independent Apoptosis Probe

Live Cell Dye

SCT238**Pack Size 1 mg****Store at -20 °C****FOR RESEARCH USE ONLY****Not for use in diagnostic procedures. Not for Human or Animal Consumption.**

Background

The BioTracker™ Apo-15 Calcium-Independent Apoptosis Probe is a calcium-independent, “turn-on” fluorescence probe emitting in the green spectrum for imaging and detection of apoptotic live cells.

Apoptosis is dysregulated in numerous diseases, from inflammatory conditions to cancer. Detection of apoptosis is a key output in assessing drug efficacy, yet the low levels of free calcium in diseased tissues limit use of annexins as apoptosis detection reagents, favoring calcium-independent fluorogenic detection. Apo-15 binds negatively charged phospholipids exposed on apoptotic cell membranes without interfering with cellular function. Apo-15 is non-fluorescent in viable cells but highly fluorescent in apoptotic cells, enabling highly specific detection of early apoptosis. Apoptotic imaging and detection with Apo-15 is broadly applicable to a wide range of cell types and experimental conditions.

Spectral Properties

Fluorescence images obtained by $\lambda_{\text{ex}} = 500 \text{ nm}$ and emission at 520 nm.

Quality Control Testing

Purity $\geq 98\%$ confirmed by HPLC

Identification confirmed by HNMR, LC-MS, and elemental analysis

Molar Mass: 1311.33 g/mol

Storage and Handling

Store BioTracker™ Apo-15 Calcium-Independent Apoptosis Probe at -20°C, desiccated and protected from light.

Note: Centrifuge vial briefly to collect contents at bottom of vial before opening.

Presentation

Lyophilized. Orange solid.

Representative Data

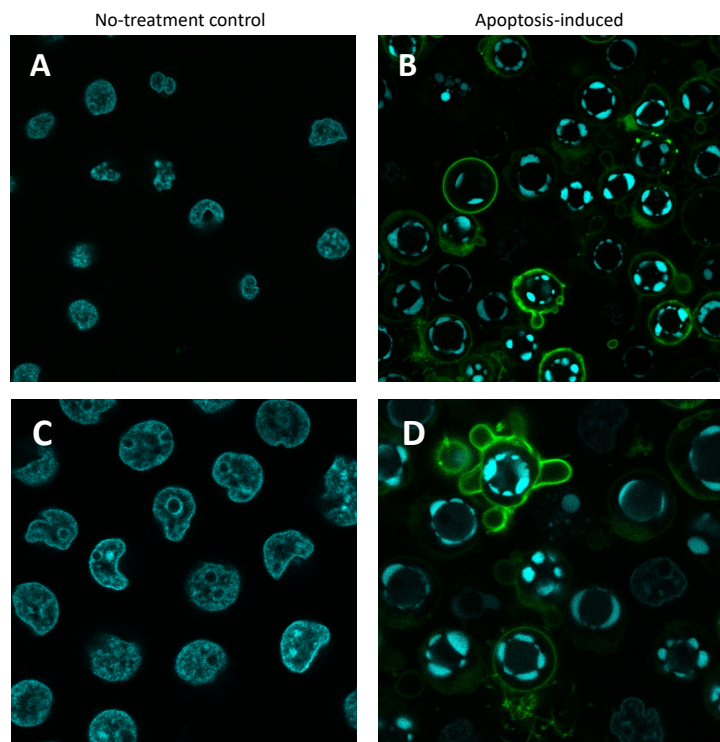


Figure 1: Confocal microscopy images of Apo-15 staining. HL-60 human leukemia cells were cultured with or without the apoptosis-inducer staurosporine (1 μ M; Cat. No. 19-123) and stained with 100 nM Apo-15 dye solution (green) and co-stained with 7 μ M Hoechst nuclear dye (cyan). Merged images show no staining in control (A, C) and staining of apoptotic cell membranes in treated samples (B, D).

Protocols

Preparing BioTracker™ Apo-15 live cell dye stock solution

1. Before opening the vial, spin down the solid to the bottom by a microcentrifuge or by a desktop centrifuge.
2. Warm the vial to room temperature. Prepare the Apo-15 probe dye stock solution by dissolving the contents of one vial (1 mg) in 763 μ L of DMSO to create a 1 mM solution.
3. Aliquot and store stock solution at -20° C or below for longer storage.

Labeling cells

1. Culture cells in an appropriate medium and vessel for fluorescence microscopy or flow cytometry
2. Prepare the Apo-15 staining solution by diluting the Apo-15 stock solution 1:10,000 to 1:1,000 in culture medium.
3. Treat cells with appropriate apoptosis inducer (example, 1 μ M staurosporine for 3 hours)
4. Remove the cell culture medium from the cells.
5. Add sufficient staining solution to cover the cells.
6. Incubate for 10-20 minutes, protected from light.
7. Observe the cells under fluorescence microscope: λ_{ex} = 500 nm, λ_{em} = 510-530 nm or detect by flow cytometry with blue/cyan (488 nm) laser.

Note: Optimal concentration must be determined by end user.

References

1. Barth N et al. A fluorogenic cyclic peptide for imaging and quantification of drug-induced apoptosis. Nat. Commun. 2020, 11:4027.

Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

Technical Assistance

Visit the tech service page at SigmaAldrich.com/techservice.

Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at SigmaAldrich.com/terms.

Contact Information

For the location of the office nearest you, go to SigmaAldrich.com/offices.

The life science business of Merck operates
as MilliporeSigma in the U.S. and Canada.

Merck, BioTracker and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2023-2024 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

Document Template 20306518 Ver 6.0

20714848 Ver 3.0, Rev 31May2024, CJ

