

Rapid and Efficient Esterification Process for FAME Profiling of Vegetable Fixed Oils and Fats

Raj Mahat, Lance Kailey, MilliporeSigma (Laramie, WY, USA)

Introduction

Vegetable fixed oils and fats have multiple valuable applications in food, cosmetics, pharmaceutical preparations, bio-fuel (biodiesel) production, and other areas.^{1, 2} In pharmaceutical industry, vegetable oils are used as an excipient to achieve improved stability and bio-availability and thus better consumer acceptability. They are used as a green solvent in drug formulations and extraction of bio-active compounds. Type I hydrogenated oils, which have high melting point and low iodine values, are used as lubricant in tablet and capsule formulation and as an alternative to hard waxes in topical pharmaceutical formulations.³ Furthermore, vegetable oil generated nanostructured lipid particles can also be used as a carrier in lipid-based formulations. Besides having nutraceutical values, vegetable oils also have a number of pharmacological and therapeutic applications, such as antimicrobial, antitumor, anti-inflammatory, and antioxidant agents.⁴

Vegetable oils and fats are liquids or semisolids which primarily consist of triacylglycerols (TAGs, ~95%), diacylglycerols (DAGs, < 5%), tocopherols and other vitamins (<1%), and phytosterols (~1%).⁵ The main component, TAGs, are made up of three fatty acyl chains substituted to a glycerol backbone. The fatty acyl chains can have 6-24 carbon atoms and be either saturated and unsaturated. Saturated fatty acids do not have a double bond while monounsaturated fatty acids have one double bond and polyunsaturated fatty acids have two or more (up to 6) double bonds.⁶ A TAG can have three of the same or different fatty acyl chains. In **Figure 1** is a typical TAG consisting of stearyl (C18:0), stearyl (C18:0), and oleyl (C18:1n9cis) fatty acyl chains shown.

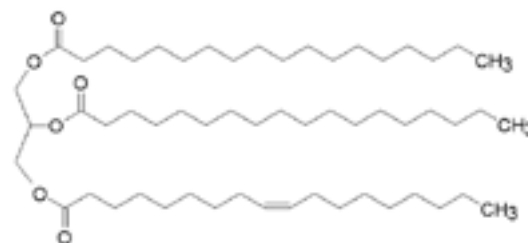


Figure 1. Structure of a typical TAG with stearyl, stearyl, and oleyl (SSO) fatty acyl chains.

Fatty acid profiling is one of the most common ways of characterizing a vegetable oil. In pharmaceutical industries, USP general chapter <401> for Fats and Fixed Oils is a common way to conduct the esterification process.⁷ However, this method suffers from a number of limitations. First, it is a two-step esterification process that uses toxic chemicals, such as boron trifluoride (BF₃). Second, and perhaps the most important one, from a practical standpoint, the method uses refluxing such that parallel sample preparation requires investment in the glasswares, such as condenser, round-bottom (RB) flasks, etc., and lab space.⁷ As a result, the method might be low-throughput and it may not be applicable for routine QC analysis.

The objective of this application note thus is to develop and validate a simple, safe, and complete alternative esterification method that is highly accurate, consistent, and capable of parallel processing for esterification of multiple vegetable oil and fat samples and also has applicability in quality control (QC) process.

Experimental

Method

Development of an alternative esterification method

A single-step, acid-catalyzed esterification process that utilizes anhydrous acetyl chloride solution (2%, v/v) in methanol to make methanolic hydrochloride reagent was improved and validated for vegetable oils.⁸

Preparation of methanolic hydrochloride reagent

To prepare the methanolic hydrochloride reagent solution, 2 mL of acetyl chloride (**Cat. No. 00990**) was mixed with 18 mL of methanol (**Cat. No. 34860**). The resulting mixture produces approximately 2 M methanolic hydrochloride (anhydrous).

Caution should be taken when adding acetyl chloride to methanol. Sudden contact with water and alcohols should be avoided. This is a highly exothermic process and can build heat and pressure if not handled

appropriately. To avoid the hazard, about 5 mL of methanol was added to a 40 mL amber vial with a PTFE septum lined cap. Acetyl chloride was added slowly with constant mixing and cooling of the vial in cold water. The pressure was vented slowly, and the vial was cooled throughout the addition of acetyl chloride. Once all the acetyl chloride was added, the rest of the methanol (~13 mL) was added into the mixture. The vial was sealed and stored free from exposure to water. The reagent can be saved for future use also if stored in a refrigerator in a properly sealed air-tight container.

Esterification reaction

About 25 mg of neat corn oil sample (**Cat. No. PHR2897**) and 2 mL of the methanolic hydrochloride reagent were mixed in a clear glass reaction vial with conical bottom. The reaction mixture was then heated in a digital two-block heater at 80 °C for 20 min. The reaction vials were then allowed to cool to room temperature. An acid-catalyzed general esterification reaction of a TAG to produce fatty acid methyl esters is shown in **Figure 2**.

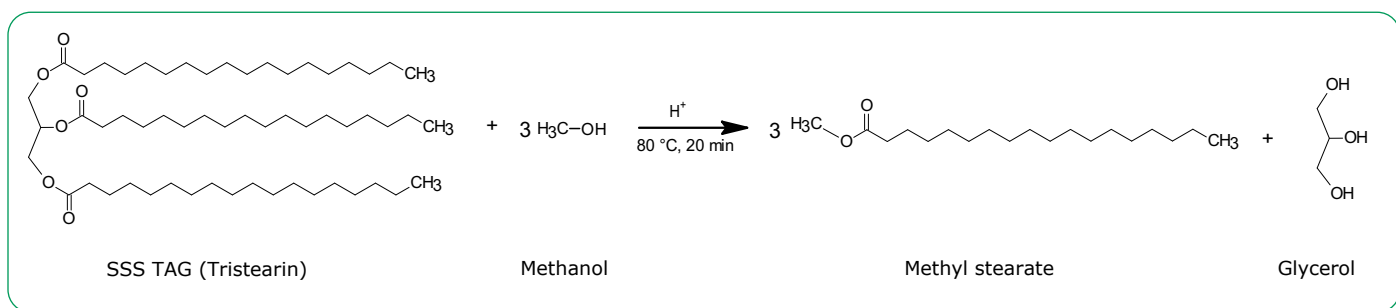


Figure 2. Acid-catalyzed methyl esterification of a TAG with stearyl (C18:0), stearyl (C18:0), and stearyl (C18:0) [SSS] fatty acyl chains to produce methyl stearate.

Extraction

To carry out the liquid-liquid extraction, 2 mL of 6% (w/v) Na₂CO₃ (**Cat. No. 222321**) solution in water was added to the cooled reaction mixture and vigorously shaken. One volume of 2 mL n-heptane (**Cat. No. 650536**) was added as the extraction solvent. After adding heptane and shaking, the mixture was allowed to stand at room temperature until the phases had completely separated. The hydrophobic fatty acid methyl esters partitioned into upper heptane phase, of which 1 mL was carefully transferred to an autosampler vial and 1 µL was injected on to a GC-FID (Agilent, 7890 A GC) equipped with a Supelco® SUPELCOWAX™ 10 column 30 m x 0.53 mm ID, 0.5 µm (**Cat. No. 25325**) for chromatographic separation. The general workflow for the esterification and detection process is shown in **Figure 3**. The details of the GC-FID detection method are presented in **Table 1** and a representative chromatogram of a corn oil FAME profile is shown in **Figure 4**.

Table 1. Chromatographic conditions used to separate FAMES from vegetable oils

Column:	SUPELCO WAX™ 10, 30 m x 0.53 mm I.D., 0.50 µm film thickness (25325)
Detector:	FID @ 280 °C
Inlet:	220° C
Column temp:	70 °C (Hold 2 min) @ 5 °C/min to 240 °C (Hold 8 min)
Flow:	4.8 mL/min
Carrier gas:	Hydrogen
Injection:	1 µL - Split injection with 20:1 split ratio
Liner:	78.5 mm x 6.3 mm, 4.0 mm I.D., Split/ Splitless, Single Taper FocusLiner™, wool packed (2879825-U)
Sample diluent:	Heptane

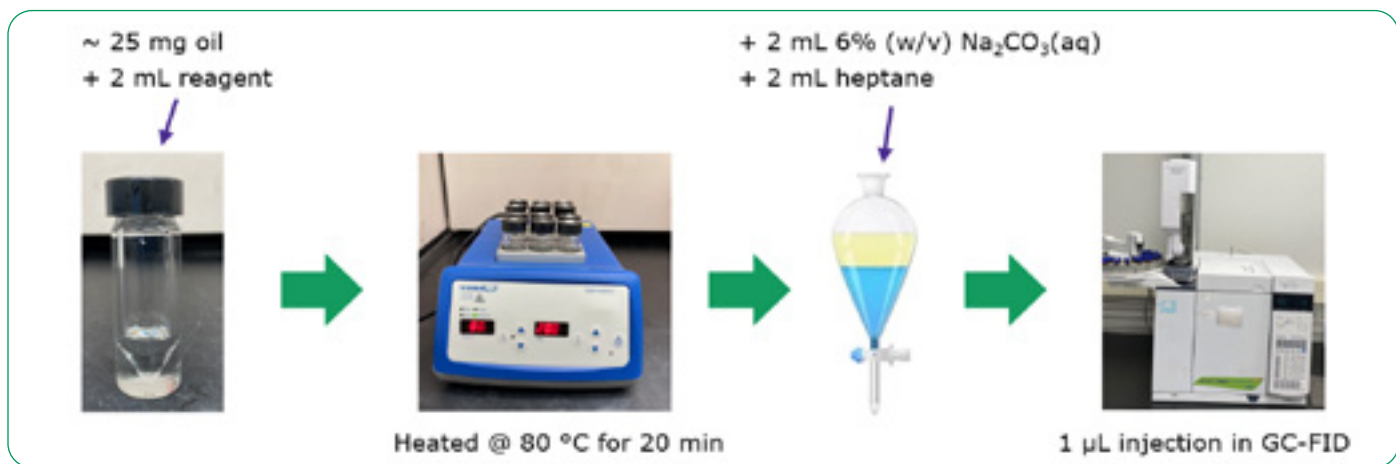


Figure 3. Representation of esterification process workflow followed by a GC-FID detection.

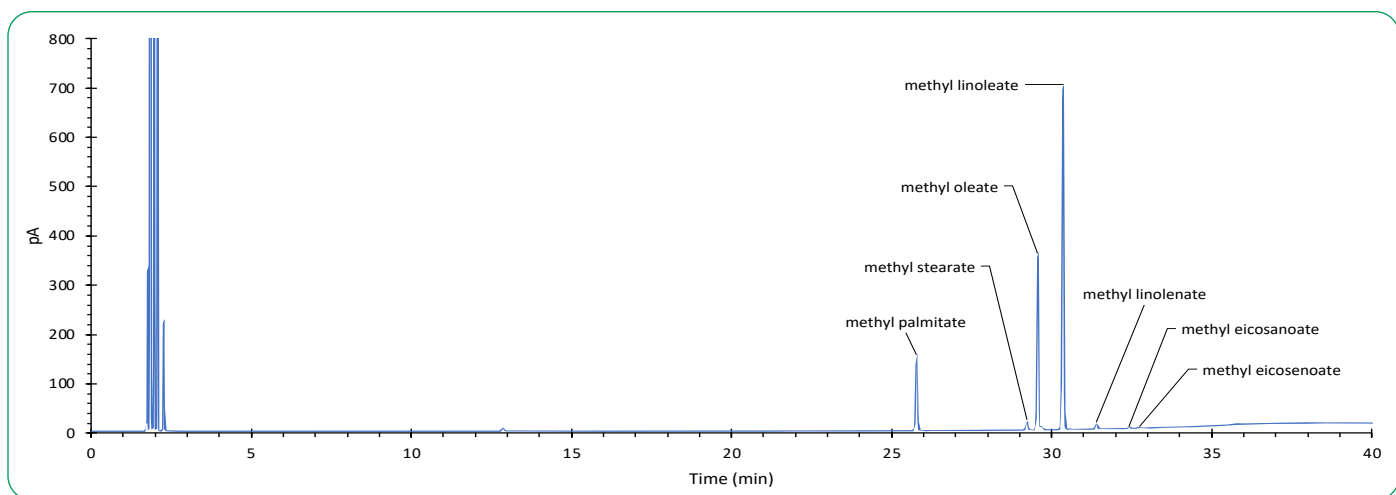


Figure 4. A representative GC-FID chromatogram showing FAME components in corn oil.

Results & Discussion

Method Validation: Quantitative esterification process

A corn oil secondary pharma standard (Cat. No. PHR2897) in five different amounts in the range of 10-35 mg was esterified and analyzed by GC-FID. The corresponding FID signals (peak area) were plotted against the amount of oil. A separately esterified oil sample having weight about 25 mg was injected 10 times to calculate precision. About 25 mg of a pharmacopeia primary standard for corn oil was esterified using the same alternative method and used as a reference standard for estimation of accuracy. Acceptance criteria for method validation parameters were referenced from USP <401> and general practices adopted while following compendial methods. Accuracy was chosen to be within 98-102%. Precision or repeatability, in term of RSD%, was set below 5% for the individual FAMES and below 1% for the ratio of methyl palmitate and methyl stearate (C16:0/C18:0).⁷ Linearity criterion was set as regression coefficient, R^2 , above 0.99. Chromatographic resolution of at least 1.5 between methyl stearate (C18:0) and methyl oleate (C18:1) peaks was the criterion for specificity.

In brief, accuracies for two predominant FAMES, methyl oleate (C18:1) and methyl linoleate (C18:2), were 101.3% and 99.8% respectively (Table 2). RSDs for all the individual FAMES were below 5% and the RSD for methyl palmitate to methyl stearate ratio (C16:0/C18:0) was 0.2% (Table 3). Linear detector responses were observed for oil samples in the amount of 10-35 mg with $R^2 > 0.99$ for all individual FAMES present. For instance, R^2 of 0.996 was found for methyl palmitate (C16:0) (Figure 5). Limit of detection (LOD) and limit of quantitation (LOQ) were estimated from LINEST function to be 2.36 mg and 7.87 mg in term of the amount of the oil sample, which correspond to 1.18 mg/mL and 3.94 mg/mL respectively in term of concentration (Table 4). The resolution between the two peaks, C18:0 and C18:1, was 3.4.

Table 2. FAME composition of corn oils (PHR2897 and pharmacopeia reference standard) follow the monograph guidelines for each individual ester. The accuracy of the esterification method was evaluated using a primary reference standard.

FAME	Shorthand notation	Supelco® PHR2897 Comp%	Primary standard Comp%	Accuracy %	Monograph Guidelines %
Methyl palmitate	C16:0	12.12	12.42		8.6-16.5
Methyl stearate	C18:0	1.60	1.63		1.0-3.3
Methyl oleate	C18:1	28.82	28.45	101.30	20-42.2
Methyl linoleate	C18:2	55.78	55.89	99.80	39.4-62
Methyl linolenate	C18:3	1.06	0.94		0.5-1.5
Methyl eicosanoate	C20:0	0.32	0.35		<0.8
Methyl eicosenoate	C20:1	0.31	0.32		<0.5

Table 3. Repeatability of the esterification method evaluated by calculating RSD% for C16:0, C18:0, and their ratio.

Repeatability estimated as RSD% for C16:0/C18:0 peak areas			
FAME	Methyl palmitate	Methyl stearate	C16:0/C18:0
Rep 1	638.635	84.665	7.54
Rep 2	636.733	84.805	7.51
Rep 3	641.531	85.060	7.54
Rep 4	632.594	83.814	7.55
Rep 1	638.635	84.665	7.54
Rep 2	636.733	84.805	7.51
Rep 3	641.531	85.060	7.54
Rep 4	632.594	83.814	7.55
Rep 5	632.928	83.917	7.54
Rep 6	643.537	85.326	7.54
Rep 7	643.548	85.137	7.56
Rep 8	639.390	84.568	7.56
Rep 9	644.542	85.393	7.55
Rep 10	641.041	84.841	7.56
Average	639.45	84.75	7.54
SD	4.27	0.54	0.01
RSD%	0.67	0.64	0.20

Efficiency of esterification reaction: Evaluation of completeness of the reaction


The completeness of the esterification process, and efficiency of the reaction, was evaluated qualitatively. An incomplete/inefficient esterification process partially hydrolyzes/esterifies TAGs producing free fatty acids and MAGs and DAGs. So, presence of such lipid classes were taken as markers for the incomplete esterification process.

For this, about 35 mg of corn oil samples were esterified with 1 and 2 mL of the methanolic hydrochloride reagent in parallel. The extractions were carried out with 2 mL of heptane at acidic pH with prior addition of 2 mL of water instead of Na₂CO₃ solution. About 500 µL of each extract was subjected to another derivatization step with an addition of 100 µL of MSTFA reagent (Cat. No. 69479) and heating for 15 min at 70 °C. This derivatization step replaces the acidic protons from the carboxylate and hydroxyl functional groups with a trimethylsilyl (TMS) moiety converting unesterified free fatty acids into -TMS ester and MAGs and DAGs into -(n)TMS ether derivatives. Here, n equals to number of hydroxyl functional group(s) in the molecular structure. A general

Table 4. Linearity, LOD, and LOQ estimation

Amount of oil (mg)	Peak Area	
9.90	236.43	
16.20	373.21	
20.80	505.27	
31.00	705.67	
34.70	770.58	
LINEST Function		
Slope	21.680	29.998
SD	0.830	20.192
R ²	0.996	17.062
LOD	2.36 mg	
LOQ	7.87 mg	

Figure 5. Calibration curve showing linear detector (GC-FID) response for the sample amount in the range of 10-35 mg of corn oil using methyl palmitate (C16:0) as the representative FAME.

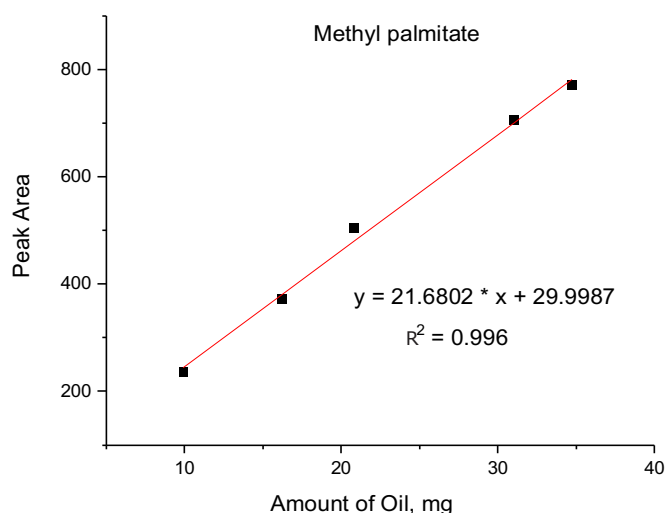

scheme of a silylation reaction for a partially hydrolyzed monoacylglyceride, such as 1-palmitin, and a free fatty acid, such as palmitic acid, is shown in Figure 6. The reaction mixture was analyzed on a single quadrupole GCMS (Agilent, 5975C MSD) in full-scan mode in the m/z range of 40-550 with a solvent delay of 3.5 min. The chromatographic separation was carried out on a SLB®-5ms (30 m x 0.25 mm x 0.25 µm, Cat. No. 28471-U) capillary column. An aliquot of 1.0 µL of the sample was injected without removal of the excess MSTFA reagent in the injector heated at 220 °C with split ratio of 10:1. Injection of silylation reagent is safe on GC columns that do not have free hydroxyl group. Said that, columns containing PEG stationary phase, such as Omegawax, Carbowax, etc., should be avoided. The GC oven was set at 70 °C for 2 min and then ramped at the rate of 5 °C/min to final temperature of 240 °C which was held for 5 min. Helium was used as carrier gas at the constant flow rate of 1 mL/min. The ionization energy of 70 eV was used to collect EI spectra at the sampling rate of 4 scans/sec. The transfer line,

ionization source, and the quadrupole analyzer were set at 240 °C, 230 °C, and 150 °C, respectively.

As shown in **Figure 7**, careful inspection of the chromatogram (Top) for the oil sample esterified with 1 mL of the methanolic hydrochloride showed presence of 1-monopalmitin-2TMS, 1-monoolein-2TMS, and 1-monolinolein-2TMS derivatives. This signified incomplete esterification process with 1 mL of the

esterification reagent. On the other hand, absence of such peaks on the bottom chromatogram suggested completeness of the process with 2 mL of the reagent. This concluded, 2 mL of the methanolic hydrochloride reagent is sufficient for complete esterification of the oil samples up to 35 mg.

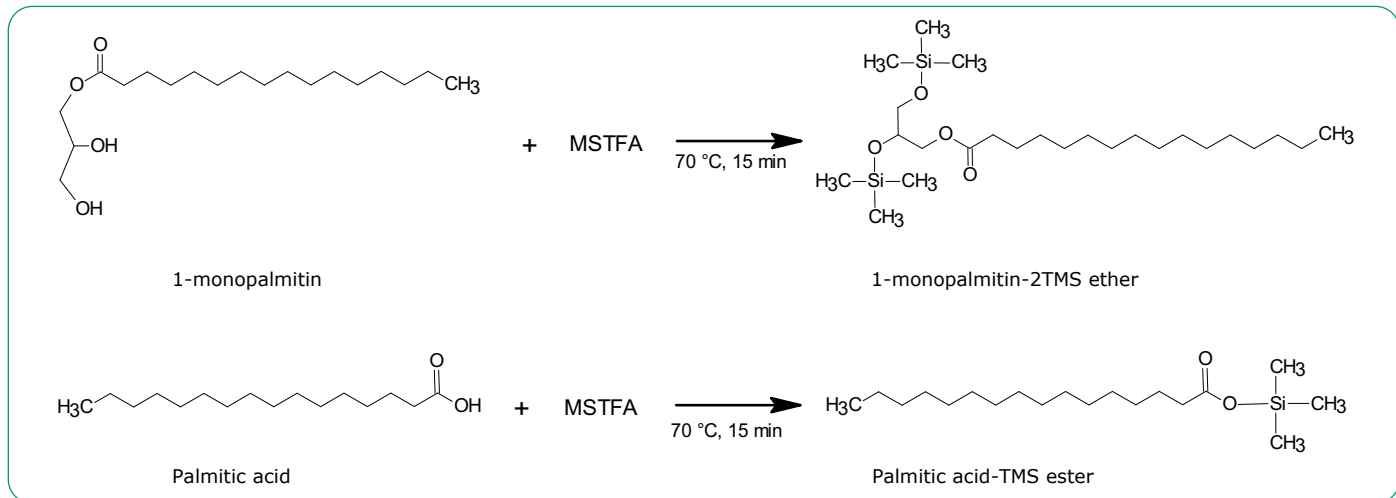


Figure 6. Trimethylsilylation derivatization of a partially esterified monoacylglycerol, such as 1-monopalmitin, to produce 1-monopalmitin-2TMS ether derivative (Top reaction). A non-esterified free fatty acid, such as palmitic acid, produces palmitic acid-TMS ester (Bottom reaction).


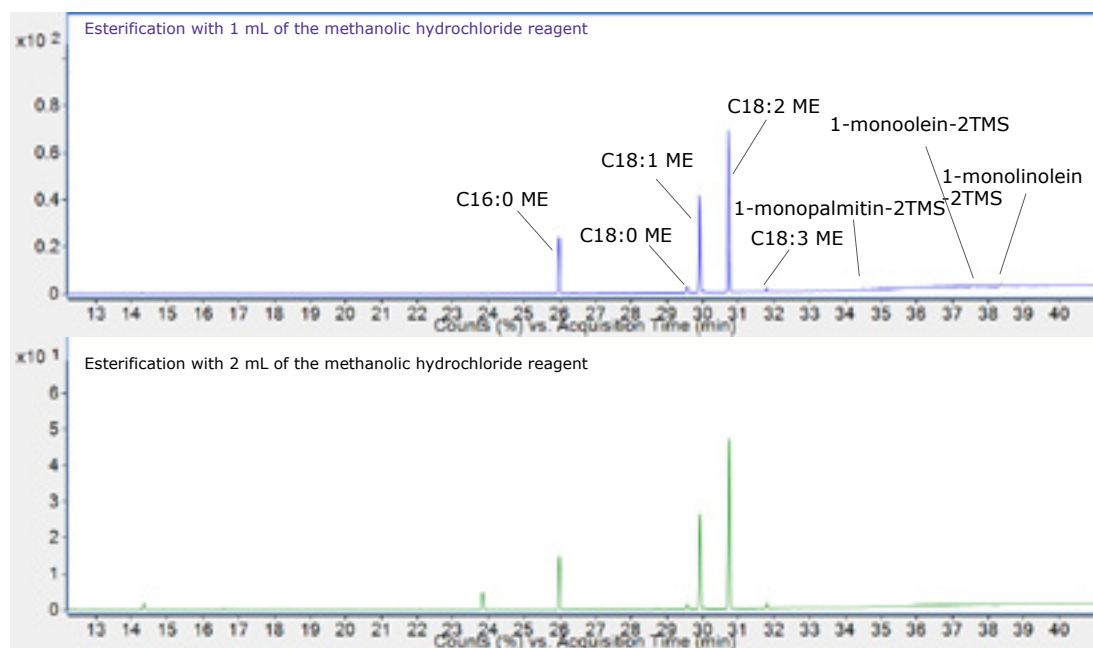


Figure 7. Chromatograms showing efficiency of the esterification process with 1 mL and 2 mL reagent.

Replication of the esterification process for analysis of olive oil samples

The esterification process was replicated for analysis of other oil samples. **Figure 8** shows the complete qualitative agreement between olive oil pharmaceutical secondary standard (**Cat. No. PHR2902**) and the pharmacopeia reference standard.

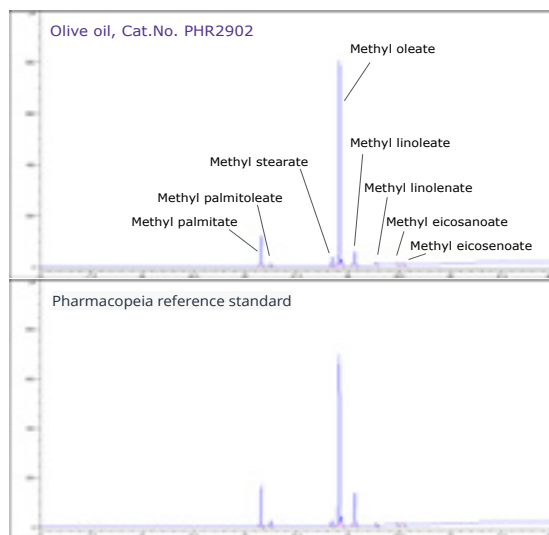

Application of the method in esterification of other vegetable oil and fat samples

The method was used in the methyl esterification of 18 fixed oil and fat samples. Replicate sample processing was carried out for each oil sample. The applicability of the method for QC of the multiple oil samples was examined by conducting the esterification of the corresponding pharmacopeia reference standards. High qualitative and quantitative FAME profile agreement between the PHR and pharmacopeia reference standards shows method's QC applicability. Summary of FAME profiles of the vegetable fixed oil and fat samples is presented in **Table 5**.

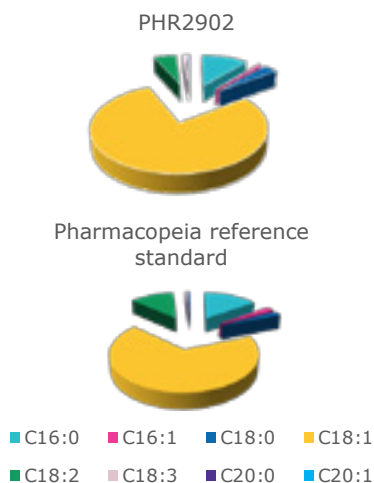


Figure 8. Comparative gas chromatograms (left) showing equivalent FAME profiles of olive oil pharmaceutical secondary standard (**Cat. No. PHR2902**) and a pharmacopeia reference standard. The pie-charts show actual FAME contents of the both standards following USP guidelines.

Table 5. A summary of FAME profiles of eighteen vegetable fixed oil and fat samples determined by esterification process that used methanolic hydrochloride reagent made from 2% (v/v) acetyl chloride solution in methanol. All the samples showed equivalent FAME profile with the corresponding pharmacopeia reference standards.

No	Cat. No.	Oil	Lot Number	Individual FAME content%							
				≤C10:0	C12:0	C14:0	C16:0	C16:1	C18:0	C18:1	C18:2
1	PHR2989	Sunflower Oil	LRAC7633				6.25		2.92	33.05	57.78
2	PHR2969	Almond Oil	LRAC7630				6.34	0.48	1.41	67.99	23.78
3	PHR2984	Palm Oil	LRAC7555		1.02	1.35	43.99		4.03	40.27	9.34
4	PHR2985	Peanut Oil	LRAC7553				9.96		2.70	55.34	24.52
5	PHR2897	Corn Oil	LRAC6778				12.12		1.60	28.82	55.77
6	PHR2902	Olive Oil	LRAC6773				10.80	0.78	3.29	78.57	5.20
7	PHR2983	Menhaden Fish Oil	LRAC7663			8.47	16.34	12.42	2.83	4.98	
8	PHR3520	Coconut Oil	LRAC6922	14.12	48.27	18.73	8.79		2.55	6.09	1.45
9	PHR2981	Hydrogenated Castor Oil	LRAC8230				1.13		9.64		
10	PHR2971	Canola Oil	LRAC7839				4.08	0.22	1.48	63.00	19.86
11	PHR2980	Flax Oil	LRAC7836				5.53		3.26	16.71	15.81
12	PHR2976	Cottonseed Oil	LRAC7835				21.24		2.61	20.42	55.73
13	PHR2987	Safflower Oil	LRAC7692				6.91		2.55	9.70	80.21
14	PHR2974	Cod Liver Oil	LRAC8228			4.77	10.00	10.06	1.97	24.07	1.78
15	PHR2970	Borage Seed Oil	LRAC8254				10.66		4.81	20.79	44.65
16	PHR2979	Fish Oil	LRAC8231			9.27	17.14	14.24	2.92	5.86	1.69
17	PHR2882	Soybean Oil	LRAD0590				11.25		3.55	19.30	56.8
18	47115-U	Lard Oil	LRAC9635			0.73	34.49	6.10	7.99	35.87	13.86

*Methyl 12-hydroxystearate is the main component (85.56%) in hydrogenated castor oil (**Cat.No. PHR2981**)

Conclusion

The developed and validated fatty acid methyl esterification method that utilizes methanolic hydrochloride reagent prepared from 2% (v/v) acetyl chloride solution in methanol is quantitative, simple, complete, and high-throughput allowing parallel sample processing. As many as 24 samples can be processed in parallel, within 20 min. The method is able to esterify various oil and fat samples including complex oil samples, such as fish oil. The method

has high QC applicability too. Strong qualitative and quantitative agreements in the results of analyses of multiple vegetable fixed oils and fats in parallel with the respective pharmacopeia reference standard confirm method's applicability in QC laboratories. This method can be used as an alternative to USP <401> in characterization/FAME profiling of vegetable oils and fats in pharmaceutical and food industry laboratories.

References

1. Enferadi Kerenkan A, Béland F, Do T-O. Chemically catalyzed oxidative cleavage of unsaturated fatty acids and their derivatives into valuable products for industrial applications: a review and perspective. *Catalysis science & technology*;6(4):971–987. <https://pubs.rsc.org/en/content/articlelanding/2016/cy/c5cy01118c>. doi:10.1021/acsomega.1c04692
2. Roy K, Poompiew N, Pongwisuthiruchte A, Potiyaraj P. Application of different vegetable oils as processing aids in industrial rubber composites: A sustainable approach. *ACS omega*. 2021;6(47):31384–31389. <http://dx.doi.org/10.1021/acsomega.1c04692>.
3. Cheikhoussef N, Cheikhoussef A. Vegetable oils as green solvents in the pharmaceutical industry. In: *Green Sustainable Process for Chemical and Environmental Engineering and Science*. Elsevier; 2021. p. 1–11. <https://doi.org/10.1016/B978-0-12-821885-3.00017-7>
4. Fagionato Masiero J, Barbosa EJ, de Oliveira Macedo L, de Souza A, Nishitani Yukuyama M, Arantes GJ, Bou-Chacra NA. Vegetable oils in pharmaceutical and cosmetic lipid-based nanocarriers preparations. *Industrial crops and products*. 2021;170(113838):113838. <http://dx.doi.org/10.1016/j.indcrop.2021.113838>
5. Hammond, E.W.; *Encyclopedia of Food Sciences and Nutrition*. 2nd edition, 2003, <https://www.sciencedirect.com/referencework/9780122270550/encyclopedia-of-food-sciences-and-nutrition>
6. Kumar A, Sharma A, C. Upadhyaya K. Vegetable oil: Nutritional and industrial perspective. *Current genomics*. 2016;17(3):230–240. <https://doi.org/10.2174/138920291766616020220107>
7. USP, USP-NF, <401> Fats and Fixed Oils, https://doi.org/10.31003/USPNF_M99160_04_01
8. Mahat R, Seebart C, Basile F, Ward NL. Global and targeted lipid analysis of *Gemmata obscuriglobus* reveals the presence of lipopolysaccharide, a signature of the classical Gram-negative outer membrane. *Journal of bacteriology*. 2016;198(2):221–236. <http://dx.doi.org/10.1128/jb.00517-15>.

Individual FAME content%									Class of FAME content (%)			
C18:3	C20:0	C20:1	C20:5	C22:0	C22:1	C22:6	C24:0	Others	Total	Total saturated	Total monounsaturated	Total polyunsaturated
									100	9.17	33.05	57.78
									100	7.75	68.47	23.78
									100	50.39	40.27	9.34
0.19	1.22	1.58		2.91	1.58				100	16.79	58.50	24.71
1.06	0.32	0.31							100	14.04	29.13	56.83
0.70	0.41	0.25							100	14.50	79.60	5.90
			14.49			12.88		27.59	100	30.32	21.53	48.15
									100	78.34	6.09	1.45
	0.41							88.82*	100			
10.09	0.40	0.87							100	5.96	64.09	29.95
58.69									100	8.79	16.71	74.50
									100	23.85	20.42	55.73
								0.63	100	9.46	9.70	80.21
0.88		12.8	9.37		7.10	10.24		6.96	100	17.10	54.03	28.87
13.98	0.33	2.39		0.29	1.33			0.77	100	16.09	24.51	58.63
			15.36			11.38		22.14	100	30.63	24.37	45.00
8.30	0.25	0.21		0.23			0.11		100	15.39	19.51	65.10
0.75								0.21	100	43.21	42.18	14.61

Featured Products

Description	Cat. No.
GC & Reference Materials	
SUPELCO WAX™ 10 Capillary GC Column, 30 m x 0.32 mm, df 0.5 µm	25325
SLB®-5ms Capillary GC Column, 30 m x 0.25 mm, df 0.25 µm	28471-U
GC Liner, Split/Splitless Type, Straight FocusLiner™ Design (wool packed)	2879825-U
Corn oil, pharmaceutical secondary standard, CRM	PHR2897
Olive oil, pharmaceutical secondary standard, CRM	PHR2902
Reagents & Accessories	
Acetyl chloride, puriss. p.a., ≥99.0%	00990
Heptane, HPLC Plus, for HPLC, GC, and residue analysis, 99%	650536
Methanol, HPLC grade, ≥99.9%	34860
MSTFA, for GC derivatization, LiChropur™, ≥98.5%	69479
Sodium carbonate, ACS reagent, anhydrous, ≥99.9%	222321

Description	Cat. No.
Oil Standards, pharmaceutical secondary standards, CRMs*	
Sunflower Oil	PHR2989
Almond Oil	PHR2969
Palm Oil	PHR2984
Peanut Oil	PHR2985
Corn Oil	PHR2897
Olive Oil	PHR2902
Menhaden Fish Oil	PHR2983
Coconut Oil	PHR3520
Hydrogenated Castor Oil	PHR2981
Canola Oil	PHR2971
Flax Oil	PHR2980
Cottonseed Oil	PHR2976
Safflower Oil	PHR2987
Cod Liver Oil	PHR2974
Borage Seed Oil	PHR2970
Fish Oil	PHR2979
Soybean Oil	PHR2882
Lard Oil (analytical standard)	47115-U

*if not indicated otherwise

See more information of Pharmaceutical Analysis & QC applications at SigmaAldrich.com/PharmaQC

Find more Analytical Reference Materials for Pharmaceutical QC at

SigmaAldrich.com/pharmastandards

To place an order or receive technical assistance

Order/Customer Service: SigmaAldrich.com/order

Technical Service: SigmaAldrich.com/techservice

SigmaAldrich.com

We have built a unique collection of life science brands with unrivalled experience in supporting your scientific advancements.

Millipore® Sigma-Aldrich® Supelco® Milli-Q® SAFC® BioReliance®

Merck KGaA
Frankfurter Strasse 250
64293 Darmstadt, Germany

