




SIGMA QUALITY CONTROL TEST PROCEDURE

Product Information

Enzymatic Assay of D-LACTIC DEHYDROGENASE (EC 1.1.1.28)

Sigma Prod. Nos. L-2011, L-2395, L-3888, and L-9636

PRINCIPLE:

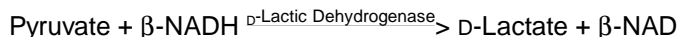

Abbreviations used:

β -NADH = β -Nicotinamide Adenine Dinucleotide, Reduced Form

β -NAD = β -Nicotinamide Adenine Dinucleotide, Oxidized Form

CONDITIONS: T = 25°C, pH = 7.0, A_{340nm}, Light path = 1 cm

METHOD: Continuous Spectrophotometric Rate Determination

REAGENTS:

- A. 100 mM Potassium Phosphate Buffer, pH 7.0 at 25°C
(Prepare 200 ml in deionized water using Potassium Phosphate, Monobasic, Anhydrous, Sigma Prod. No. P-5379. Adjust to pH 7.0 at 25°C with 1 M KOH.)
- B. 11 mM β -Nicotinamide Adenine Dinucleotide, Reduced Form Solution (β -NADH)
(Prepare 1 ml in cold deionized water using β -Nicotinamide Adenine Dinucleotide, Reduced Form, Disodium Salt, Sigma Prod. No. N-8129. **PREPARE FRESH.**)
- C. 20 mM Sodium Pyruvate Solution (Pyruvate)
(Prepare 1.0 ml in cold deionized water using Pyruvic Acid, Sodium Salt, Sigma Prod. No. P-2256.)
- D. 1.0% (w/v) Bovine Serum Albumin Solution (BSA)
(Prepare 50 ml in Reagent A using Albumin, Bovine, Sigma Prod. No. A-4503 or equivalent.)
- E. D-Lactic Dehydrogenase Enzyme Solution
(Immediately before use, prepare a solution containing 0.3 - 0.60 unit/ml of D-Lactic Dehydrogenase in cold Reagent D. **PREPARE FRESH.**)

**Enzymatic Assay of D-LACTIC DEHYDROGENASE
(EC 1.1.1.28)
Sigma Prod. Nos. L-2011, L-2395, L-3888, and L-9636**

PROCEDURE:

Pipette (in milliliters) the following reagents into suitable cuvettes:

	<u>Test</u>	<u>Blank</u>
Reagent A (Buffer)	2.50	2.50
Reagent B (β-NADH)	0.05	0.05
Reagent C (Pyruvate)	0.10	0.10

Mix by inversion and equilibrate to 25°C. Monitor the $A_{340\text{nm}}$ until constant, using a suitably thermostatted spectrophotometer. Then add:

Reagent D (BSA)	-----	0.05
Reagent E (Enzyme Solution)	0.05	-----

Immediately mix by inversion and record the decrease in $A_{340\text{nm}}$ for approximately 5 minutes. Obtain the $\Delta A_{340\text{nm}}/\text{minute}$ using the maximum linear rate for both the Test and Blank.

CALCULATIONS:

$$\text{Units/ml enzyme} = \frac{(\Delta A_{340\text{nm}}/\text{min Test} - \Delta A_{340\text{nm}}/\text{min Blank})(2.7)(\text{df})}{(6.22)(0.05)}$$

2.7 = Total volume (in milliliters) of the assay

df = Dilution

6.22 = Millimolar extinction coefficient of β-NADH at 340 nm

0.05 = Volume (in milliliters) of assay

$$\text{Units/mg solid} = \frac{\text{units/ml enzyme}}{\text{mg solid/ml enzyme}}$$

$$\text{Units/mg protein} = \frac{\text{units/ml enzyme}}{\text{mg protein/ml enzyme}}$$

UNIT DEFINITION:

One unit will reduce 1.0 μmole of pyruvate to D-lactate per minute at pH 7.0 at 25°C.

FINAL ASSAY CONCENTRATION:

In a 2.75 ml reaction mix, the final concentrations are 94 mM potassium phosphate, 0.20 mM β-nicotinamide adenine dinucleotide, reduced form, 0.74 mM pyruvate, 0.015 - 0.03 unit D-lactic dehydrogenase.

**Enzymatic Assay of D-LACTIC DEHYDROGENASE
(EC 1.1.1.28)
Sigma Prod. Nos. L-2011, L-2395, L-3888, and L-9636**

NOTES:

1. This assay is based on the cited reference.
2. Where Sigma Product or Stock numbers are specified, equivalent reagents may be substituted.

Sigma warrants that the above procedure information is currently utilized at Sigma and that all Sigma-Aldrich, Inc. products conform to the information in this and other Sigma-Aldrich, Inc. publications. Purchaser must determine the suitability of the information and product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packaging slip.