EnglishMillex®-SV Filter Unit (for low-binding applications)Cat. No. SLSV025LS (50/box)

- Single use only
- Sterile
- Non-pyrogenic

Introduction

The bidirectionally supported Millex®-SV device is a low-binding, syringe-operated filter unit for the filtration of aqueous, proteinaceous solutions where minimal sample loss is desired. It will remove particles, precipitates, and undissolved powders larger than 5.0 microns (µm). The sterile Millex®-SV filter unit is non-pyrogenic and non-toxic. This single-use device consists of a 5.0 µm Durapore® membrane filter sealed in a plastic housing. Typical applications include particle or aggregate removal from protein solutions such as monoclonal antibodies, extracted cellular or tissue-based solutions, hormones, and drugs.

Chemical Compatibility


The Millex®-SV filter unit is compatible with most aqueous solutions. It may be used to filter the agents listed below. This guide has been developed from technical publications, materials suppliers, and laboratory tests, and is believed to be reliable. However, because of variability in temperature, concentrations, duration of exposure, and other factors outside of our control which may affect the use of the unit, no warranty is given or is to be implied with respect to such information. Agents not listed below should be tested with the Millex®-SV filter unit prior to use.

Acetic acid (glacial)	Butyl alcohol	Hydrofluoric acid	Paraldehyde
Acetic acid (5%)	Carbon tetrachloride	Hydrogen (gas)	Pentane
Aliphatic ethers	Ethyl alcohol	Hydrogen peroxide (3%)	Petroleum based oils
Ammonium hydroxide (6 N)	Ethylene glycol	Hypo (photo)	Petroleum ether
Amyl acetate	Formaldehyde	Isobutyl alcohol	Phenol (0.5%)
Amyl alcohol	Freon® solvent, TF or PCA	Isopropyl alcohol	Pyridine
Benzyl alcohol (1%)	Glycerine (glycerol)	Methyl alcohol	Silicone oils
Boric acid	Helium	Mineral spirits	
Brine (sea water)	Hexane	Nitrogen	

Directions for Use

WARNINGS:

- Do to ensure sterility, do not use this product if the package is damaged.
- Do not use this product as an in-line filter for intravenous administration; it was not designed for long-term continuous use.
- Do not use with syringes smaller than 10 mL because pressures in excess of the maximum pressure rating may be reached, potentially causing damage to the filter unit and/or personal injury.
- CAUTIONS:
 - Do not resterilize or reuse this filter unit, as Merck Millipore Ltd. cannot assure the sterility, integrity, and performance beyond a single use.
 - Do not use the Millex®-SV unit at temperatures above 45 °C (113 °F).
 - Do not use the same Millex®-SV unit to filter solutions in both directions.
 - Do not use the Millex®-SV unit to filter emulsions or suspensions.
 - Do not use the Millex®-SV unit to filter 5 mg or less of active drug materials unless binding studies have been performed.

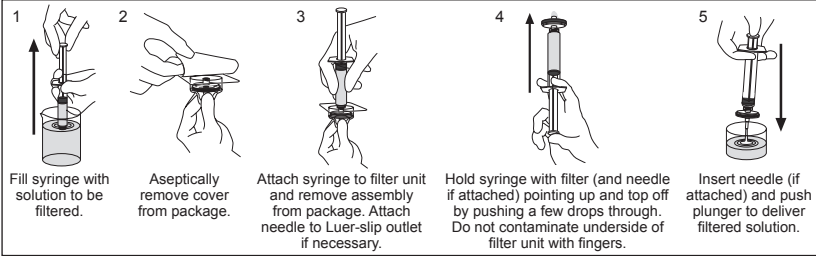Specifications

Membrane	Hydrophilic Durapore® PVDF (polyvinylidene fluoride), type SV, pore size 5.0 µm
	PVC (polyvinyl chloride)
Housing	
Inlet to outlet	25 mm (0.98 in.)
Diameter	29 mm (1.14 in.)
Filtration area	4 cm² (0.62 in²)
Temperature limit	45 °C (113 °F)
Pressure limit at 25 °C	5.2 bar (75 psi) inlet and differential
Filtration volume	1–100 mL
Hold-up volume	≤ 0.1 mL after air purge
Sterilization method	Ethylene oxide gas
Connections	Female Luer-Lok™ inlet, male Luer-slip outlet
Flow rate at 2.1 bar (30 psi), 25 °C	≥ 600 mL/min

Notice

The information in this document is subject to change without notice and should not be construed as a commitment by Merck Millipore Ltd. ("Millipore") or an affiliate. Neither Merck Millipore Ltd. nor any of its affiliates assumes responsibility for any errors that may appear in this document.

Technical Assistance

For more information, contact the office nearest you. In the U.S., call 1-800-MILLIPORE (1-800-645-5476). Outside the U.S., go to our web site at www.millipore.com/offices for up-to-date worldwide contact information. You can also visit the tech service page on our web site at www.millipore.com/techservice.

Standard Product Warranty

The applicable warranty for the products listed in this publication may be found at www.millipore.com/terms (within the "Terms and Conditions of Sale" applicable to your purchase transaction).

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