

Data Sheet

# BioTracker™ MultiTASQ G-quadruplex G4 Cell Probe

**SCT247****Pack Size: 1 mg****Store at -20 °C****FOR RESEARCH USE ONLY****Not for use in diagnostic procedures. Not for human or animal consumption.**

## Background

G-quadruplexes are four-stranded structures comprising stacked G tetrads (G4) formed within certain guanine-rich nucleic acid sequences. G-quadruplexes are enriched in open chromatin regions but are also present in the nucleolus, ribosomal regions, and cytoplasm. The formation of G4 structures is dynamic and cell-type specific, suggestive of interaction with transcription factors and other proteins.<sup>1</sup> For this reason G-quadruplexes have garnered intense interest for their potential functions in transcriptional regulation.

The BioTracker™ MultiTASQ G-quadruplex G4 Cell Probe is optimized for cellular detection of G-quadruplexes. MultiTASQ was designed utilizing the template-assembled synthetic G-quartets (TASQ) concept as biomimetic ligand for G-quadruplexes, demonstrating high specificity for G4. MultiTASQ is water-soluble and exhibits a high degree of intracellular accessibility. MultiTASQ detects ribosomal RNA-G4 associated with perinuclear rough ER regions to a greater degree compared to BG4 antibody. MultiTASQ is an alkynylated probe that allows for background-free detection via click chemistry and is suitable for G4 capture chemi-precipitation and click imaging applications. MultiTASQ is compatible with live-cell incubation. The BioTracker™ MultiTASQ G4 Cell Probe is a uniquely designed, versatile tool for G4 detection and analysis.

## Source

BioTracker™ MultiTASQ G-quadruplex G4 Cell Probe (SCT247) does not contain genetically modified organisms.

## Spectral Properties

This probe is non-fluorescent. Probe may be detected using click (copper-catalyzed cycloaddition) imaging with appropriate click chemistry reagents.

## Quality Control Testing

Purity: ≥ 98% confirmed by HPLC, HNMR, LC-MS and elemental analysis


Molar Mass: 1800 g/mol

## Storage and Handling



Store BioTracker™ MultiTASQ G-quadruplex G4 Cell Probe at -20 °C, desiccated and protected from light.

**Note:** Centrifuge vial briefly to collect contents at bottom of vial before opening.


## Representative Data

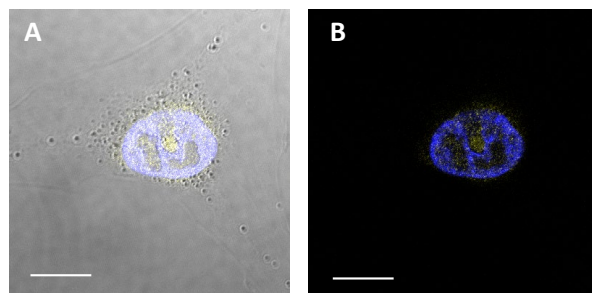


**Figure 1.** Confocal microscopy images of G4 foci in HeLa cells detected with anti-DNA G-quadruplex structure BG4 antibody (MABE917) and anti-FLAG antibody (MAB3118, yellow), co-stained with DAPI nuclear dye (blue). (**A, B**) No BG4 control vs (**C, D**) BG4 stained cells. (**A, C**) overlay with brightfield images. White arrows indicate G4 foci in nucleus. Scale bars: 10  $\mu$ M

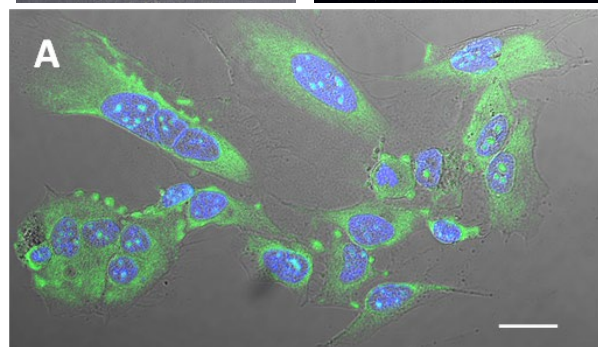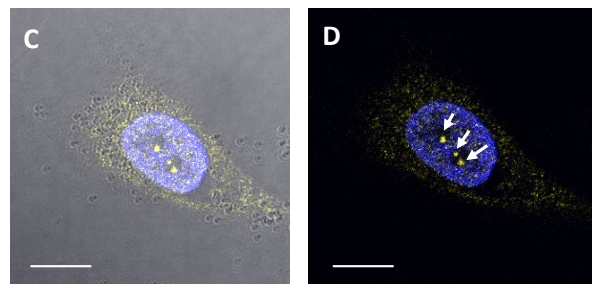

**Figure 2.** Confocal microscopy images of G-quadruplexes detected with MultiTASQ. HeLa cells were incubated with MultiTASQ then fixed and incubated with Atto-488-conjugated azide (72709, green) and co-stained with DAPI nuclear dye (blue). (**A**) overlay with brightfield; (**B**) overlay. Nucleolar, perinuclear, and cytoplasmic foci and diffuse staining is observed. Scale bars: 25  $\mu$ M

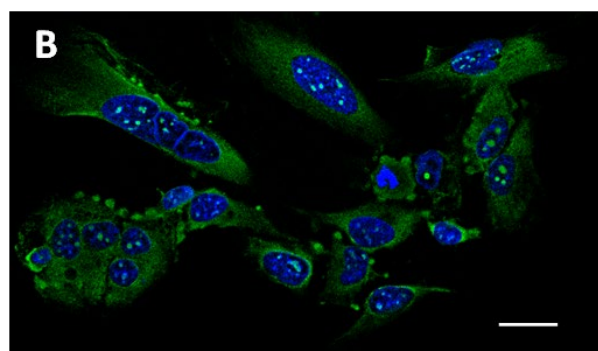

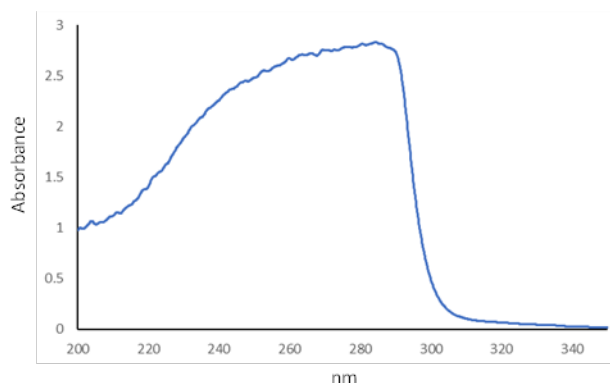

**Figure 3.** Absorption spectrum of BioTracker™ MultiTASQ G-quadruplex G4 Cell Probe (SCT247) in water.

## Protocols

### Preparing BioTracker™ MultiTASQ G4 Cell Probe stock solution

1. Before opening the vial, spin down the solid to the bottom by a microcentrifuge or by a desktop centrifuge.
2. Warm the vial to the room temperature. Prepare the MultiTASQ (Molecular Weight: 1800 g/mol) probe stock solution by dissolving the contents of one vial (1 mg) in 1.8 mL of water to create a 1 mM solution.
3. Aliquot and store stock solution at  $-20^{\circ}\text{C}$  or below for longer-term storage.

### Preparing Click Reaction Buffer

Prepare a solution of PBS containing 0.05% (v/v) IGEPAL® CA-630 (I3021), 10 mM sodium ascorbate (A4034), and 1 mM copper sulfate (209198).

### Labeling cells

1. Culture cells in an appropriate medium and vessel for fluorescence microscopy.
2. Prepare the MultiTASQ staining solution by diluting the MultiTASQ stock solution 1:1000 in growth medium (1  $\mu\text{M}$  final concentration).
3. Add sufficient staining solution to cover the cells and incubate for 24 hours at  $37^{\circ}\text{C}$  in 5%  $\text{CO}_2$  humidified incubator.
4. Wash cells 3 times for 5 minutes with PBS.
5. Fix cells (for example, ice-cold methanol for 10 minutes).
6. Wash cells 3 times for 5 minutes with PBS.
7. Incubate cells with 1  $\mu\text{M}$  fluorophore-conjugated azide (for example, Atto488-azide, 72709) in click reaction buffer for 1 hour at ambient temperature.
8. Wash cells for 5 minutes with PBS. If desired, counterstain with DAPI (1  $\mu\text{g}/\text{mL}$ ) for 10 minutes at ambient temperature.
9. Image cells on fluorescence microscope at wavelengths appropriate for fluorophores used.

## Presentation

Lyophilized. White solid.

## References

1. Nat. Chem. 2021, 13(7): 626-633.
2. ACS Chem Biol. 2021, 16(5): 905-914.

---

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

## Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://SigmaAldrich.com/techservice).

## Terms and Conditions of Sale

Warranty, use restrictions, and other conditions of sale may be found at [SigmaAldrich.com/terms](https://SigmaAldrich.com/terms).

## Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://SigmaAldrich.com/offices).

The life science business of Merck operates  
as MilliporeSigma in the U.S. and Canada.

Merck, BioTracker and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2008-2024 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

Document Template 20306518 Ver 6.0

23087738 Ver 1.0, Rev 14AUG2024, LB

